PRODUCT INFORMATION & MANUAL

Human/Monkey IL-5 Valukine™ ELISA

Catalog Number: VAL125

For the quantitative determination of natural and recombinant
human/monkey IL-5 concentrations

For research use only.
Not for diagnostic or therapeutic procedures.

Bio-Techne China Co. Ltd
P: +86 (21) 52380373 P: 8009881270 F: +86 (21) 52381001
info.cn@bio-techne.com

Please refer to the kit label for expiry date.
Novus kits are guaranteed for 3 months from date of receipt

Version 202403.5

TABLE OF CONTENTS

I. BACKGROUND	2
II. OVERVIEW	3
III. ADVANTAGES	4
IV. EXPERIMENT	7
V. KIT COMPONENTS AND STORAGE	8
VI. PREPARATION	10
VII. ASSAY PROCEDURE	12
VIII. REFERENCES	14

I. BACKGROUND

Interleukin-5 (IL-5) is a secreted glycoprotein that belongs to the alpha-helical group of cytokines (1-3). Unlike other family members, it is present as a covalently linked antiparallel dimer (4, 5). The cDNA for human IL-5 encodes a signal peptide and a 115 amino acid (aa) mature protein. Mature human IL-5 shares 70%, 70%, 62%, 71%, 70% and 66%, aa sequence identity with mouse, rat, canine, equine, feline and porcine IL-5, respectively and shows cross-reactivity with mouse IL-5. IL-5 is primarily produced by CD4⁺ Th2 cells, but also by activated eosinophils, mast cells, EBV-transformed B cells, Reed-Sternberg cells in Hodgkin's disease, and IL-2-stimulated invariant natural killer T cells (iNKT) (1-3, 6-8). IL-5 increases production and mobilization of eosinophils and CD34⁺ progenitors from the bone marrow and causes maturation of eosinophil precursors outside the bone marrow (1, 6, 9, 10).

The receptor for human IL-5, mainly expressed by eosinophils, but also found on basophils and mast cells, consists of a unique ligand-binding subunit (IL-5 R alpha) and a shared signal-transducing subunit, beta c (3, 6, 11). IL-5 R alpha first binds IL-5 at low affinity, then associates with preformed beta c dimers, forming a high-affinity receptor (12). IL-5 also binds proteoglycans, potentially enhancing its activity (13). Soluble forms of IL-5 R alpha antagonize IL-5 and can be found in vivo (10, 14).

In humans, IL-5 primarily affects cells of the eosinophilic lineage, and promotes their differentiation, maturation, activation, migration and survival, while in mice IL-5 also enhances Ig class switching and release from B1 cells (1-3, 9, 10, 15, 16). IL-5 also promotes differentiation of basophils and primes them for histamine and leukotriene release (17).

II. OVERVIEW

A. PRINCIPLE OF THE ASSAY

This assay employs the quantitative sandwich enzyme immunoassay technique. An antibody specific for IL-5 has been pre-coated onto a microplate. Standards and samples are pipetted into the wells and any IL-5 present is bound by the immobilized antibody. After washing away any unbound substances, a biotinylated detection antibody specific for IL-5 is pipetted into the wells. After washing away any unbound substances, Streptavidin-HRP is pipetted into the wells. Following a wash to remove any unbound reagent, TMB substrate solution (Chromogenic agent) is added to the wells and color develops in proportion to the amount of IL-5 bound in the initial step. The color development is stopped, and the intensity of the color is measured.

B. LIMITATIONS OF THE PROCEDURE

- **FOR RESEARCH USE ONLY. NOT FOR USE IN DIAGNOSTIC PROCEDURES.**
- This kit is suitable for cell culture supernate and human/monkey serum.
- The kit should not be used beyond the expiration date on the kit label.
- Do not mix or substitute reagents with those from other lots or sources.
- If samples generate values higher than the highest standard, dilute the samples with Reagent Diluent (1×) and repeat the assay.
- Any variation in operator, pipetting technique, washing technique, incubation time or temperature, and kit age can cause variation in binding.

III. ADVANTAGES

A. PRECISION

Intra-assay Precision (Precision within an assay)

Three samples were tested twenty times on one plate to assess intra-assay precision.

Inter-assay Precision (Precision between assays)

Three samples were tested in twenty separate assays to assess inter-assay precision.

	Intra-assay Precision			Inter-assay Precision		
Sample	1	2	3	1	2	3
Mean (pg/mL)	8.7	34.0	124.3	8.8	34.5	125.6
Standard Deviation	80.3	1.0	4.1	0.5	1.3	7.5
CV%	3.0	2.9	3.3	5.4	3.7	6.0

B. RECOVERY

The recovery of IL-5 spiked to different levels throughout the range of the assay in cell culture media was evaluated. The recovery ranged from 88.3 to 99.1% with an average of 94.6%.

The recovery of IL-5 spiked to different levels throughout the range of the assay in human serum was evaluated. The recovery ranged from 83.0 to 89.1% with an average of 85.7%.

The recovery of IL-5 spiked to different levels throughout the range of the assay in monkey serum was evaluated. The recovery ranged from 84.1 to 90.7% with an average of 87.8%.

C. SENSITIVITY

The minimum detectable dose (MDD) of IL-5 is typically less than 1.99 pg/mL.

The MDD was determined by adding two standard deviations to the mean optical density value of twenty zero standard replicates and calculating the corresponding concentration.

D. CALIBRATION

This immunoassay is calibrated against a highly purified *sf 21*-expressed recombinant human IL-5 produced at R&D Systems®.

E. LINEARITY

To assess the linearity of the assay, different samples were containing or spiked with high concentrations of IL-5 and diluted with Reagent Diluent (1×) to produce samples with values within the dynamic range of the assay.

Dilution		Human		Monkey	
		Cell culture media (n=3)	Serum (n=3)	Cell culture media (n=3)	Serum (n=3)
1:2	Average % of Expected	99	100	101	102
	Range (%)	98 - 100	98 - 102	98 - 104	99 - 104
1:4	Average % of Expected	101	98	101	101
	Range (%)	98 - 102	96 - 99	98 - 102	99 - 104
1:8	Average % of Expected	100	96	102	96
	Range (%)	96 - 103	93 - 99	96 - 103	96 - 97
1:16	Average % of Expected	100	91	102	82
	Range (%)	96 - 103	88 - 93	96 - 108	74 - 87

F. SAMPLE VALUES

Human serum - Five human serum samples were evaluated for the presence of IL-5 in this assay. All samples measured less than the lowest IL-5 standard, 4.7 pg/mL.

Monkey serum - Three monkey serum samples were evaluated for the presence of IL-5 in this assay. All samples measured less than the lowest IL-5 standard, 4.7 pg/mL.

G. SPECIFICITY


The following factors prepared at 50 ng/mL were assayed and exhibited no cross-reactivity or interference.

Recombinant human:	Other recombinants:
GM-CSF	Bovine IL-5
IL-3	Canine IL-5
IL-5 R α	Equine IL-5
IL-5 R β	Feline IL-5
	Mouse IL-5
	Porcine IL-5
	Rabbit IL-5
	Rat IL-5

IV. EXPERIMENT

EXAMPLE STANDARD

The standard curve is provided for demonstration only. A standard curve should be generated for each set of samples assayed.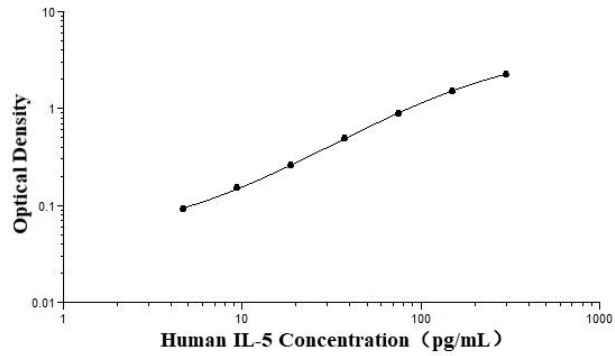

pg/mL	O.D.	Average	Corrected
0	0.036 0.043	0.040	-
4.7	0.090 0.092	0.091	0.051
9.4	0.148 0.152	0.150	0.110
18.8	0.259 0.263	0.261	0.221
37.5	0.481 0.491	0.486	0.446
75	0.893 0.891	0.892	0.852
150	1.507 1.540	1.524	1.484
300	2.241 2.272	2.257	2.217

V. KIT COMPONENTS AND STORAGE

A. MATERIALS PROVIDED

Parts	Description	Size
Human IL-5 Microplate	96 well polystyrene microplate (12 strips of 8 wells) coated with an antibody against IL-5.	1 plate
Human IL-5 Standard	Recombinant IL-5 in a buffered protein base; lyophilized. Refer to the vial label for reconstitution volume.	2 vials
Human IL-5 Detection Antibody	Biotinylated IL-5 antibody, lyophilized. Refer to the vial label for reconstitution volume.	1 vial
Streptavidin-HRP B (40×)	40× Streptavidin conjugated to horseradish peroxidase.	1 vial
Reagent Diluent Concentrate (10×)	A 10× concentrated buffered protein base used to dilute standard, samples, detection antibody and HRP.	1 vial
Wash Buffer Concentrate (25×)	A 25× concentrated solution of buffered surfactant with preservatives.	1 vial
TMB Substrate	TMB ELISA Substrate Solution/ TMB Substrate Solution	1 vial
Stop Solution	2 N sulfuric acid.	1 vial
Plate Sealers	Adhesive strips.	3 strips

B. STORAGE

Unopened Kit	Store at 2-8 °C. Do not use past kit expiration date.	
Opened/ Reconstituted Reagents	Streptavidin-HRP B	May be stored for up to 1 month at 2-8 °C.*
	Wash Buffer (1×)	
	TMB Substrate	
	Stop Solution	
	Standard	Prepare fresh for each assay. Standards may be stored for up to 1 month at -20°C*.
	Detection Antibody	Aliquot and store for up to 1 month at -20 °C in a manual defrost freezer.*
	Reagent Diluent Concentrate (10×)	May be stored for up to 1 month at 2-8 °C.* Use and discard diluted Reagent Diluent (1×). Prepare fresh for each assay.
Microplate Wells	Return unused wells to the foil pouch containing the desiccant pack, reseal along entire edge of zip-seal. May be stored for up to 1 month at 2-8 °C.*	

* Provided this is within the expiration date of the kit.

C. OTHER SUPPLIES REQUIRED

- ◆ Microplate reader capable of measuring absorbance at 450 nm, with the correction wavelength set at 540 nm or 570 nm.
- ◆ Pipettes and pipette tips.
- ◆ Deionized or distilled water.
- ◆ Squirt bottle, manifold dispenser, or automated microplate washer.
- ◆ Test tubes for dilution of standards.
- ◆ 100 mL and 500 mL graduated cylinder.

D. PRECAUTION

- ◆ Some components in this kit contain a preservative which may cause an allergic skin reaction. Avoid breathing mist.
- ◆ The Stop Solution provided with this kit is an acid solution. Wear protective gloves, clothing, eye, and face protection. Wash hands thoroughly after handling.

VI. PREPARATION

A. SAMPLE COLLECTION AND STORAGE

Cell Culture Supernates - Remove particulates by centrifugation and assay immediately or aliquot and store samples at ≤ -20 °C. Avoid repeated freeze-thaw cycles. Samples may require dilution with Reagent Diluent (1×).

Serum - Use a serum separator tube (SST) and allow samples to clot for 30 minutes at room temperature before centrifugation for 15 minutes at 1000 × g. Remove serum and assay immediately or aliquot and store samples at ≤ -20 °C. Avoid repeated freeze-thaw cycles. Samples may require dilution with Reagent Diluent (1×).

B. SAMPLE PREPARATION

Monkey serum samples recommend a 2-fold dilution prior to the assay. A suggested 2-fold dilution is 100 μ L of sample + 100 μ L of Reagent Diluent (1×). Optimal dilutions should be determined by the end user.

C. REAGENT PREPARATION

Note: *Bring all reagents to room temperature before use.*

Wash Buffer (1×) - If crystals have formed in the concentrate, warm to room temperature and mix gently until the crystals have completely dissolved. Dilute 20 mL of Wash Buffer Concentrate (25×) into deionized or distilled water to prepare 500 mL of Wash Buffer (1×).

Reagent Diluent (1×) - Use deionized or distilled water to prepare Reagent Diluent (1×).

Detection Antibody (1×) - **Centrifuge briefly before opening. Reconstitution volume refer to vial label to prepare Detection Antibody (100×).** Allow the Detection Antibody to sit for a minimum of 15 minutes with gentle agitation prior to making dilutions. Aliquot and store if needed. Dilute to Detection Antibody (1×) with Reagent Diluent (1×). Prepare at least 15 minutes prior to use.


Streptavidin-HRP B (1×) - **Centrifuge briefly before opening.** Dilute to the working concentration specified on the vial label using Reagent Diluent (1×).

IL-5 Standard - **Centrifuge briefly before opening. Refer to the vial label for the reconstitution volume* using Reagent Diluent (1×).** This reconstitution produces a

stock solution of 3000 pg/mL. Allow the standard to sit for a minimum of 15 minutes with gentle agitation prior to making dilutions.

*If you have any question, please seek help from our Technical Support.

Pipette 900 μ L of Reagent Diluent (1 \times) into the 300 pg/mL tube. Pipette 500 μ L of the appropriate Reagent Diluent (1 \times) into the remaining tubes. Use the stock solution to produce a dilution series (below). Mix each tube thoroughly before the next transfer. The 300 pg/mL standard serves as the high standard. The Reagent Diluent (1 \times) serves as the zero standard (0 pg/mL).

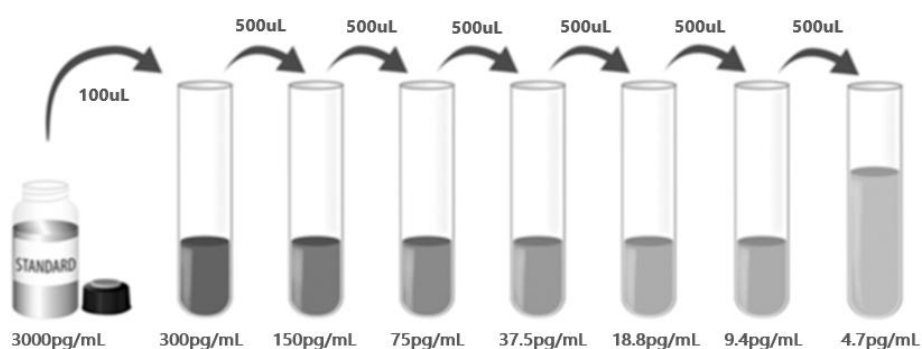D. TECHNICAL HINTS

- When mixing or reconstituting protein solutions, always avoid foaming.
- To avoid cross-contamination, change pipette tips between additions of each standard level, between sample additions, and between reagent additions. Also, use separate reservoirs for each reagent.
- It is recommended that the samples be pipetted within 15 minutes.
- To ensure accurate results, proper adhesion of plate sealers during incubation steps is necessary.
- TMB Substrate should remain colorless until added to the plate. Keep TMB Substrate protected from light. TMB Substrate should change from colorless to gradations of blue.
- Stop Solution should be added to the plate in the same order as the TMB substrate. The color developed in the wells will turn from blue to yellow upon addition of the Stop Solution. Wells that are green in color indicate that the Stop Solution has not mixed thoroughly with the TMB substrate.

VII. ASSAY PROCEDURE

Note: Bring all reagents and samples to room temperature before use. It is recommended that all samples and standards be assayed in duplicate.

1. Prepare all reagents and working standards as directed in the previous sections.
2. Remove excess microplate strips from the plate frame, return them to the foil pouch containing the desiccant pack, and reseal.
3. Add 100 μ L of standard or sample per well. Cover with the adhesive strip provided. **Incubate for 2 hours at room temperature.** A plate layout is provided for a record of standards and samples assayed.
4. Aspirate each well and wash, repeating the process three times for a total of four washes. Wash by filling each well with Wash Buffer (400 μ L) using a squirt bottle, manifold dispenser, or autowasher. Complete removal of liquid at each step is essential to good performance. After the last wash, remove any remaining Wash Buffer by aspirating or decanting. Invert the plate and blot it against clean paper towels.
5. Add 100 μ L of the Detection Antibody (1 \times) diluted in Reagent Diluent (1 \times), to each well. Cover with a new adhesive strip and **incubate 2 hours at room temperature.**
6. Repeat the aspiration/wash as in step 4.
7. Add 100 μ L of the working dilution of Streptavidin-HRP B to each well. Cover the plate and **incubate for 20 minutes at room temperature. Protect from light.**
8. Repeat the aspiration/wash as in step 4.
9. Add 100 μ L of TMB Substrate to each well. **Incubate for 20 minutes at room temperature. Protect from light.**
10. Add 50 μ L of Stop Solution to each well. Gently tap the plate to ensure thorough mixing.
11. Determine the optical density of each well within 10 minutes, using a microplate reader set to 450 nm. If wavelength correction is available, set to 540 nm or 570 nm. If wavelength correction is not available, subtract readings at 540 nm or 570 nm from the readings at 450 nm. This subtraction will correct for optical imperfections in the plate. Readings made directly at 450 nm without correction may be higher and less accurate.
12. **CALCULATION OF RESULTS**

Average the duplicate readings for each standard and sample and subtract the average zero standard optical density (O.D.). Create a standard curve by reducing the data using computer software capable of generating a four parameter logistic (4-PL) curve-fit. As an alternative, construct a standard curve by plotting the mean absorbance for each standard on the y-axis against the concentration on the x-axis and draw a best fit curve through the points on the graph. The data may be linearized by plotting the log of the IL-5 concentrations versus the log of the O.D. and the best fit line can be determined by regression analysis. This procedure will produce an adequate but less precise fit of the data. If samples have been diluted, the concentration read from the standard curve must be multiplied by the dilution factor.

VIII. REFERENCES

1. Rosenberg, H. F. et al. (2007) *J. Allergy Clin. Immunol.* 119:1303.
2. Elsas, P.X. and M. I. G. Elsas (2007) *Curr. Med. Chem.* 14:1925.
3. Martinez-Moczygamba, M. and D. P. Huston (2003) *J. Allergy Clin. Immunol.* 112:653.
4. Minamitake, Y. et al. (1990) *J. Biochem.* 107:292.
5. McKenzie, A. N. et al. (1991) *Mol. Immunol.* 28:155.
6. Shakoory, B. et al. (2004) *J. Interferon Cytokine Res.* 24:271.
7. Lalani, T. et al. (1999) *Ann. Allergy Asthma Immunol.* 82:317.
8. Sakuishi, K. et al. (2007) *J. Immunol.* 179:3452.
9. Clutterbuck, E. J. et al. (1989) *Blood* 73:1504.
10. Cameron, L. et al. (2000) *J. Immunol.* 164:1538.
11. Tavernier, J. et al. (1991) *Cell* 66:1175.
12. Zaks-Zilberman, M. et al. (2008) *J. Biol. Chem.* 283:13398.
13. Lipscombe, R. et al. (1998) *J. Leukocyte Biol.* 63:342.
14. Tavernier, J. et al. (2000) *Blood* 95:1600.
15. Kopf, M. et al. (1996) *Immunity* 4:15.
16. Horikawa, K. and K. Takatsu (2006) *Immunology* 118:497.
17. Denburg, J. A. et al. (1991) *Blood* 77:1462

PLATE LAYOUT

Use this plate layout to record standards and samples assayed.

1																	
2																	
3																	
4																	
5																	
6																	
7																	
8																	
9																	
10																	
11																	
12																	
	A	B	C	D	E	F	G	H									

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