

ELISpot

Human* IL-17

Catalog Number EL317

For the quantitative determination of the frequency of cells releasing human IL-17.

*This kit also recognizes primate IL-17.

This package insert must be read in its entirety before using this product.
For research use only. Not for use in diagnostic procedures.

TABLE OF CONTENTS

SECTION	PAGE
INTRODUCTION	1
PRINCIPLE OF THE ASSAY	2
LIMITATIONS OF THE PROCEDURE	3
TECHNICAL HINTS	3
ELISPOT SCHEMATIC	4
MATERIALS PROVIDED & STORAGE CONDITIONS	5
OTHER SUPPLIES REQUIRED	5
PRECAUTIONS	6
REAGENT PREPARATION	6
SAMPLE PREPARATION	6
ASSAY PROCEDURE	7
CALCULATION OF RESULTS	8
REPRODUCIBILITY DATA	8
SPECIFICITY	8
TROUBLESHOOTING GUIDE	9
REFERENCES	10

MANUFACTURED AND DISTRIBUTED BY:

USA & Canada | R&D Systems, Inc.

614 McKinley Place NE, Minneapolis, MN 55413, USA
TEL: (800) 343-7475 (612) 379-2956 FAX: (612) 656-4400
E-MAIL: info@RnDSystems.com

DISTRIBUTED BY:

UK & Europe | R&D Systems Europe, Ltd.

19 Barton Lane, Abingdon Science Park, Abingdon OX14 3NB, UK
TEL: +44 (0)1235 529449 FAX: +44 (0)1235 533420
E-MAIL: info@RnDSystems.co.uk

China | R&D Systems China Co., Ltd.

24A1 Hua Min Empire Plaza, 726 West Yan An Road, Shanghai PRC 200050
TEL: +86 (21) 52380373 FAX: +86 (21) 52371001
E-MAIL: info@RnDSystemsChina.com.cn

INTRODUCTION

Human Interleukin 17 (IL-17; also known as IL-17A) is a 30-35 kDa variably glycosylated homodimeric protein that belongs to a unique family of cysteine-knot related proteins (1-3). Its sequence was originally isolated from an activated rodent hybridoma and termed CTLA-8 (2-5). It is synthesized as a 155 amino acid (aa) precursor that contains a 23 aa signal sequence and a 15 kDa, 132 aa mature segment (6, 7). Although there are two intrachain disulfide bonds that create a ring reminiscent of those found in cysteine-knot proteins, the actual closed knot structure does not appear to form (1, 2). IL-17 has one potential N-linked glycosylation site. Mature human IL-17 shares 97%, 61%, and 60% aa sequence identity to rhesus macaque, mouse, and rat IL-17, respectively (4, 5, 8). It also shows limited aa sequence identity to other human IL-17 family members. In particular, it shows 34% (9), 38% (9), 34% (10) and 26% (11) aa sequence identity to IL-17B, C, D, and E, respectively, and 49% aa sequence identity to IL-17F with which it is most homologous (12). While rodent and human mature sequences show modest aa sequence identity, human IL-17 is active on both mouse and rat cells (4, 13). The cells principally known to produce IL-17 are the memory CD4⁺ T cells. In addition, CD8⁺ T cells as well as $\alpha\beta$ TCR⁺ CD4⁺CD8⁻ T cells, neutrophils (PMNs) and eosinophils have also been reported to express mRNA transcripts for IL-17 (3, 5, 14-16).

A high affinity receptor for human IL-17 (IL-17 R) has been reported and found to be ubiquitously expressed. It is synthesized as a 160 kDa, type I transmembrane glycoprotein that bears no resemblance to members of the cytokine, TNF or immunoglobulin receptor superfamily (2, 14, 17). It is synthesized as an 866 aa precursor that contains a 27 aa signal sequence, a 2931 aa extracellular region, a 21 aa transmembrane segment, and a 525 aa cytoplasmic domain (17). The extracellular region has no distinguishing aa motifs or domains. IL-17 R signalling reportedly utilizes multiple classes of MAP kinases (1), and also activates the JAK/STAT family of transcription factors (14).

Functionally, IL-17 is best known for its participation in the recruitment and survival of neutrophils (3, 14, 18). Its induction is believed to be the result of antigen stimulation of dendritic cells, resulting in IL-23 secretion. In a TCR-independent event, IL-23 can induce T cell production of IL-17 (3). Once secreted, IL-17 in the bone marrow appears to induce stromal fibroblast expression of both G-CSF and SCF (membrane form), an effect that increases PMN differentiation and production. IL-17 may complement this by directly blocking neutrophil apoptosis, promoting greater circulating PMN numbers (18). In the tissues, IL-17 appears to promote neutrophil extravasation, principally through its effects on macrophages and endothelial cells. On macrophages, IL-17 induces TNF- α , IL-1 β , and IL-6 production (19). TNF- α and IL-1 β then act on local endothelial cells to induce G-CSF secretion, an effect that is potentiated by IL-17 (20). IL-17 further contributes to PMN influx by inducing endothelial cells to release CXC chemokines and nitric oxide production, which may increase vascular permeability (3, 13).

The Human IL-17 ELISpot assay is designed for the detection of human IL-17 secreting cells at the single cell level, and it can be used to quantitate the frequency of human IL-17 secreting cells. ELISpot assays are well suited for monitoring immune responses to various stimuli, treatments and therapies, and they have been used for the quantitation of antigen-specific CD4⁺ and/or CD8⁺ T cell responses. Other methods for the assessment of antigen-specific T cell responses, such as the chromium release assays with quantitation by limiting dilution, are tedious, and require previous *in vitro* expansion of T cells for several days. These assays typically are not suitable for measuring infrequent T cell responses that occur at less than 1 in 1000. ELISpot assays are highly reproducible and sensitive, and can be used to measure responses with frequencies well below 1 in 100,000. ELISpot assays do not require prior *in vitro* expansion of T cells, and they are suitable for high-throughput analysis using only small volumes of primary cells. As such, ELISpot assays are useful tools for research in areas as diverse as antigen recognition, vaccine development, cytokine secretion and the monitoring of various clinical trials.

PRINCIPLE OF THE ASSAY

The enzyme-linked immunospot (ELISpot) assay was originally developed for the detection of individual B cells secreting antigen-specific antibodies (21, 22). This method has since been adapted for the detection of individual cells secreting specific cytokines or other antigens (23, 24). ELISpot assays employ the quantitative sandwich enzyme-linked immunosorbent assay (ELISA) technique.

A polyclonal antibody specific for human IL-17 has been pre-coated onto a polyvinylidene difluoride (PVDF)-backed microplate. Appropriately stimulated cells are pipetted into the wells and the microplate is placed into a humidified 37 °C CO₂ incubator for a specified period of time. During this incubation period, the immobilized antibody in the immediate vicinity of the secreting cells binds secreted IL-17. After washing away any cells and unbound substances, a biotinylated polyclonal antibody specific for human IL-17 is added to the wells. Following a wash to remove any unbound biotinylated antibody, alkaline-phosphatase conjugated to streptavidin is added. Unbound enzyme is subsequently removed by washing and a substrate solution (BCIP/NBT) is added. A blue-black colored precipitate forms at the sites of cytokine localization and appears as spots, with each individual spot representing an individual IL-17 secreting cell. The spots can be counted with an automated ELISpot reader system or manually using a stereomicroscope.



LIMITATIONS OF THE PROCEDURE

- FOR RESEARCH USE ONLY. NOT FOR USE IN DIAGNOSTIC PROCEDURES.
- The kit should not be used beyond the expiration date on the kit label.
- Do not mix or substitute reagents with those from other lots or sources.
- Any variation in pipetting and washing techniques, incubation time or temperature, or kit age can cause variation in the density of spots, intensity of specific staining, and background levels.


TECHNICAL HINTS

- To minimize edge effect, place the microplate (bottom down) onto a piece of aluminum foil (about 4 x 6 inches). Add cells, cover the microplate with the lid, and shape the foil around the edges of the microplate. The foil may be left on the microplate for the rest of the experimental procedure and removed after the BCIP/NBT Substrate has been washed off.
- Do not remove the flexible plastic underdrain on the bottom of the microplate before or during incubation and development. It may damage the PVDF membrane filter. The underdrain cover may be removed only after completing the incubation with the BCIP/NBT Substrate.
- Do not touch PVDF membrane filters with pipette tips when pipetting cells and reagents to avoid damage to the membrane.
- Upon completing the experiment, do not dry the microplate at a temperature above 37 °C. It may cause the PVDF membrane filters to crack.
- The 96-well microplate provided in this kit is not sterile. Due to the short incubation period and the presence of antibiotics in the culture media, microbial contamination has not been shown to be an issue with this ELISpot procedure.
- This kit is designed for single use only. The layout of the assay should be carefully planned to maximize the use of the provided microplate and reagents.
- The controls listed are recommended for each ELISpot experiment:
 - Positive Control** - Use recombinant human IL-17.
 - Unstimulated/Negative Control** - Use the same number of unstimulated cells as stimulated cells.
 - Background Control** - Use sterile culture media.
 - Detection Antibody Control** - Substitute phosphate buffered saline for detection antibody.


ELISPOT SCHEMATIC

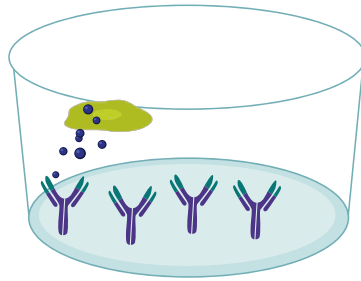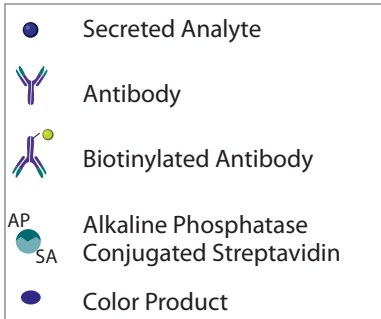


Incubate IL-17 secreting cells in an antibody-coated well.

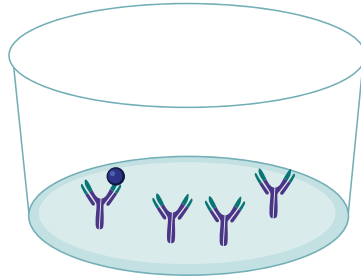

Remove cells by washing. Secreted IL-17 is captured by the immobilized antibody.

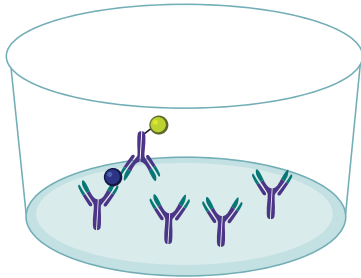


Incubate with biotinylated anti-IL-17 antibody.

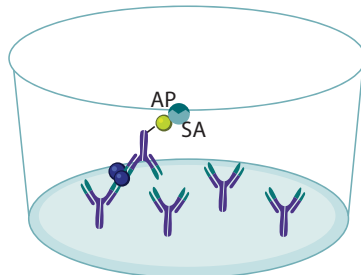

Remove unbound biotinylated antibody by washing. Incubate with alkaline phosphatase conjugated streptavidin.

Add BCIP/NBT Substrate

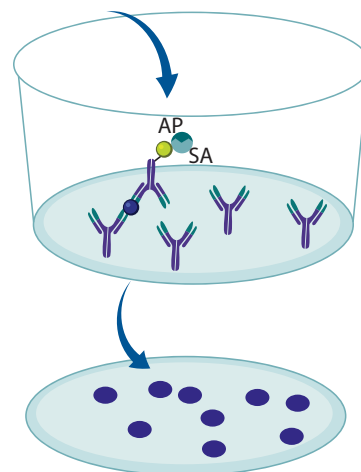

Wash to remove unbound enzyme.

Add substrate and monitor the formation of colored spots. Analyze using either an ELISpot or dissection microscope.

MATERIALS PROVIDED & STORAGE CONDITIONS

Store the unopened kit at 2-8 °C. Do not use past kit expiration date.

Note: *This kit is validated for single use only. Results obtained using previously opened or reconstituted reagents may not be reliable.*

PART	PART #	DESCRIPTION
Human IL-17 Microplate	892727	96-well PVDF-backed microplate coated with a polyclonal antibody specific for human IL-17.
Human IL-17 Detection Antibody Concentrate	892728	150 µL of a 120X concentrated solution of biotinylated polyclonal antibody specific for human IL-17 with preservatives.
Streptavidin-AP Concentrate A	895358	150 µL of a 120X concentrated solution of Streptavidin conjugated to Alkaline Phosphatase with preservatives.
Dilution Buffer 1	895307	12 mL of a buffer for diluting Human IL-17 Detection Antibody Concentrate with preservatives.
Dilution Buffer 2	895354	12 mL of a buffer for diluting Streptavidin-AP Concentrate A with preservatives.
Wash Buffer Concentrate	895308	50 mL of a 10X concentrated solution of a buffered surfactant with preservative.
BCIP/NBT Substrate	895867	12 mL of a stabilized mixture of 5-Bromo-4-Chloro-3' Indolylphosphate p-Toluidine Salt (BCIP) and Nitro Blue Tetrazolium Chloride (NBT).
Human IL-17 Positive Control	892729	3 ng of recombinant human IL-17 with preservatives; lyophilized.

OTHER SUPPLIES REQUIRED

- Dissection microscope or an ELISpot reader.
- Pipettes and pipette tips.
- Deionized water.
- Squirt bottle, manifold dispenser, or automated microplate washer.
- 500 mL graduated cylinder.
- 37 °C CO₂ incubator.
- Sterile culture media.

PRECAUTIONS

Some components of this kit contain sodium azide, which may react with lead and copper plumbing to form explosive metallic azides. Flush with large volumes of water during disposal.

BCIP/NBT is toxic if swallowed, in contact with skin, or if inhaled. It is a highly flammable liquid and vapor may cause serious irritation and damage to organs. Do not eat, drink, or smoke when using this product. Do not breathe fumes. Use only in a well-ventilated area. Keep away from heat, sparks, open flames, and hot surfaces. Keep the container tightly closed.

Some components in this kit contain a preservative which may cause an allergic skin reaction. Avoid breathing mist.

Wear protective gloves, clothing, eye, and face protection. Wash hands thoroughly after handling. Refer to the SDS on our website prior to use.

REAGENT PREPARATION

Bring all reagents to room temperature, except the Human IL-17 Detection Antibody Concentrate and Dilution Buffer 1, which should remain at 2-8 °C.

Wash Buffer - If crystals have formed in the concentrate, warm to room temperature and mix gently until the crystals have completely dissolved. To prepare Wash Buffer, add 50 mL of Wash Buffer Concentrate to 450 mL of deionized water and mix well.

Human IL-17 Positive Control - Reconstitute the lyophilized Human IL-17 Positive Control with 250 µL of culture medium that is used to incubate cells.

Detection Antibody Mixture - Tap or vortex the vial to release reagent collected in the cap. Transfer 100 µL of Human IL-17 Detection Antibody Concentrate into the vial labeled Dilution Buffer 1 and mix well. **For optimal performance, prepare the Detection Antibody immediately before use.**

Streptavidin-AP Concentrate A - Tap or vortex the vial to release reagent collected in the cap. Transfer 100 µL of Streptavidin-AP Concentrate A into the vial labeled Dilution Buffer 2 and mix well. **For optimal performance, prepare the Streptavidin-AP immediately before use.**

SAMPLE PREPARATION

The types of effector and responder cells used, method of cell separation, mode of stimulation, and length of incubation are to be determined by each investigator.

ASSAY PROCEDURE

Bring all reagents to room temperature, except the diluted Detection Antibody Mixture which should remain at 2-8 °C. All samples and controls should be assayed at least in duplicate.

1. Fill all wells in the microplate with 200 μL of sterile culture media and incubate for approximately 20 minutes at room temperature.
2. When cells are ready to be plated, aspirate the culture media from the wells. Immediately add 100 μL of the appropriate cells or controls to each well (see Technical Hints for appropriate controls).
3. Incubate cells in a humidified 37 °C CO₂ incubator. Optimal incubation time for each stimulus should be determined by the investigator. **Do not disturb the cells during the incubation period.**
4. Aspirate each well and wash, repeating the process three times for a total of four washes. Wash by filling each well with Wash Buffer (250-300 μL) using a squirt bottle, manifold dispenser, or autowasher. Complete removal of liquid at each step is essential to good performance. After the last wash, remove any remaining Wash Buffer by aspirating or decanting. Invert the plate and blot it against clean paper towels.
Note: *Adjust the height of the prongs of the manifold dispenser or autowasher to prevent damage to the membranes.*
5. Add 100 μL of Detection Antibody Mixture into each well, and incubate overnight at 2-8 °C. Alternatively, incubation with detection antibodies can be done for 2 hours at room temperature on a rocking platform.
6. Repeat the wash procedure described in step 4.
7. Add 100 μL of diluted Streptavidin-AP Concentrate A into each well, and incubate for 2 hours at room temperature.
8. Repeat the wash procedure described in step 4.
9. Add 100 μL of BCIP/NBT Substrate into each well, and incubate for 1 hour at room temperature. **Protect from light.**
10. Decant the BCIP/NBT Substrate solution from the microplate and rinse the microplate with deionized water. Invert the microplate and tap to remove excess water. Remove the flexible plastic underdrain from the bottom of the microplate, wipe the bottom of the plate thoroughly with paper towels and dry completely either at room temperature (60-90 minutes) or 37 °C (15-30 minutes).

CALCULATION OF RESULTS

The developed microplate can be analyzed by counting spots using either a dissection microscope or an ELISpot reader. Specific spots are round and have a dark center with slightly fuzzy edges. Quantitation of results can be done, for example, by calculating the number of spot forming cells (SFC) per number of cells added to the well.

REPRODUCIBILITY DATA

Human peripheral blood mononuclear cells (2.0×10^6 cells/mL) were stimulated with $4.0 \mu\text{g/mL}$ of Concanavalin A overnight at 37°C in a $5\% \text{CO}_2$ incubator. Samples were assayed in eight wells ($100 \mu\text{L/well}$) according to the procedure and analyzed with a dissection microscope.

	Number of Spots Counted
Well	Human
1	47
2	49
3	41
4	53
5	40
6	40
7	39
8	36

Rhesus and cynomolgus macaque peripheral blood mononuclear cells (2.0×10^6 cells/mL) were stimulated with $0.5 \mu\text{g/mL}$ of calcium ionomycin and $0.05 \mu\text{g/mL}$ of PMA overnight at 37°C in a $5\% \text{CO}_2$ incubator. Samples were assayed in eight wells ($100 \mu\text{L/well}$) according to the procedure and analyzed with a dissection microscope.

	Number of Spots Counted	
Well	Rhesus Macaque	Cynomolgus Macaque
1	121	202
2	127	233
3	116	232
4	117	234
5	122	196
6	117	223
7	116	235
8	127	213

SPECIFICITY

This kit detects human and primate IL-17 and does not cross-react with human IL-17B, IL-17C, or IL-17E. Low level ($\sim 5\%$) cross-reactivity has been observed with human IL-17F.

TROUBLESHOOTING GUIDE

OBSERVATION	PROBLEM	CORRECTIVE ACTION
Following the incubation with BCIP/NBT Substrate and rinsing the microplate with deionized water, the dark blue background color of the filter membrane attenuates visualization and quantitation of spots.	The membrane is wet.	Microplates cannot be analyzed accurately until the PVDF filter membranes are completely dry. Wait until the membrane becomes dry (typically 15-30 minutes at 37 °C or 60-90 minutes at room temperature).
The number of spots in the wells that contained the cells is high, but their contrast as well as intensity of staining in the Positive Control wells is low.	Underdevelopment; perhaps the result of using Streptavidin-AP and/or BCIP/NBT solutions that have not been brought to room temperature.	Warm the appropriate reagents to room temperature before adding them to the wells.
The number of spots in the wells that contained cells is lower than expected whereas Positive Control wells turned blue-black.	Cell stimulation problem.	Ensure that reagents used to stimulate the cytokine release from the cells retained their biological activity. One way to check is to perform immunocytochemistry on fixed cells after stimulation.
	Too few cells were added to the wells.	Increase the number of cells added per well.
Following incubation with BCIP/NBT Substrate and drying the microplate, the density of the spots makes them difficult to quantify.	Too many cells were added to the wells.	Make dilutions of cells (<i>i.e.</i> , 1×10^6 , 5×10^5 , 1×10^5 , 5×10^4 , 1×10^4 cells per well) to determine the optimal number of cells that will result in formation of distinct spots.

REFERENCES

1. Aggarwal, S. and A.L. Gurney (2002) *J. Leukoc. Biol.* **71**:1.
2. Moseley, T.A. *et al.* (2003) *Cytokine Growth Factor Rev.* **14**:155.
3. Kolls, J.K. and A. Linden (2004) *Immunity* **21**:467.
4. Kennedy, J. *et al.* (1996) *J. Interf. Cytokine Res.* **16**:611.
5. Rouvier, E. *et al.* (1993) *J. Immunol.* **150**:5445.
6. Yao, Z. *et al.* (1995) *J. Immunol.* **155**:5483.
7. Fossiez, F. *et al.* (1996) *J. Exp. Med.* **183**:2593.
8. GenBank Accession #: XP_001106391.
9. Li, H. *et al.* (2000) *Proc. Natl. Acad. Sci. USA* **97**:773.
10. Starnes, T. *et al.* (2002) *J. Acquir. Immune Defic. Syndr. Hum. Retrovirol.* **169**:642.
11. Lee, J. *et al.* (2001) *J. Biol. Chem.* **276**:1660.
12. Starnes, T. *et al.* (2001) *J. Immunol.* **167**:4137.
13. Miljkovic, D.J. *et al.* (2003) *Cell. Mol. Life Sci.* **60**:518.
14. Witowski, J. *et al.* (2004) *Cell. Mol. Life Sci.* **61**:567.
15. Ferretti, S. *et al.* (2003) *J. Immunol.* **170**:2106.
16. Shin, H.C. *et al.* (1998) *Cytokine* **10**:841.
17. Yao, Z. *et al.* (1997) *Cytokine* **9**:794.
18. Schwarzenberger, P. *et al.* (2000) *J. Immunol.* **164**:4783.
19. Jovanovic, D.V. *et al.* (1998) *J. Immunol.* **160**:3513.
20. Numasaki, M. *et al.* (2004) *Immunol. Lett.* **95**:97.
21. Czerkinsky, C.C. *et al.* (1983) *J. Immunol. Methods* **65**:109.
22. Sedgwick, J.D. and P.G. Holt (1983) *J. Immunol. Methods* **57**:301.
23. Czerkinsky, C.C. *et al.* (1984) *J. Immunol. Methods* **72**:489.
24. Helms, T. *et al.* (2000) *J. Immunol.* **164**:3723.

All trademarks and registered trademarks are the property of their respective owners.