

# Luminex<sup>®</sup> Performance Assay

## Human XL Cytokine Fixed Panel

Catalog Number LKTM014

For the simultaneous quantitative determination of multiple human cytokine concentrations in cell culture supernates, serum, plasma, and platelet-poor plasma.

This package insert must be read in its entirety before using this product.  
For research use only. Not for use in diagnostic procedures.

# TABLE OF CONTENTS

SECTION	PAGE
INTRODUCTION .....	1
PRINCIPLE OF THE ASSAY .....	2
LIMITATIONS OF THE PROCEDURE .....	2
MATERIALS PROVIDED & STORAGE CONDITIONS .....	3
OTHER SUPPLIES REQUIRED .....	4
PRECAUTIONS.....	4
TECHNICAL HINTS.....	4
SAMPLE COLLECTION AND STORAGE .....	5
SAMPLE PREPARATION.....	5
REAGENT PREPARATION .....	6
DILUTED MICROPARTICLE COCKTAIL PREPARATION.....	7
DILUTED BIOTIN-ANTIBODY COCKTAIL PREPARATION .....	7
STREPTAVIDIN-PE PREPARATION .....	7
INSTRUMENT SETTINGS .....	8
ASSAY PROCEDURE .....	9
ASSAY PROCEDURE SUMMARY .....	10
CALCULATION OF RESULTS .....	11
CALIBRATION .....	11
PRECISION .....	12
ACCURACY .....	13
SPECIFICITY.....	15
PLATE LAYOUT .....	17

## Manufactured and Distributed by:

### USA R&D Systems, Inc.

614 McKinley Place NE, Minneapolis, MN 55413

TEL: 800 343 7475 612 379 2956

FAX: 612 656 4400

E-MAIL: info@bio-techne.com

## Distributed by:

### Europe | Middle East | Africa Bio-Techne Ltd.

19 Barton Lane, Abingdon Science Park

Abingdon OX14 3NB, UK

TEL: +44 (0)1235 529449

FAX: +44 (0)1235 533420

E-MAIL: info.emea@bio-techne.com

### China Bio-Techne China Co., Ltd.

Unit 1901, Tower 3, Raffles City Changning Office,

1193 Changning Road, Shanghai PRC 200051

TEL: +86 (21) 52380373 (400) 821-3475

FAX: +86 (21) 52371001

E-MAIL: info.cn@bio-techne.com

## INTRODUCTION

Cytokines are intercellular signaling proteins released from a wide variety of cells and tissues. They play an integral role in regulating growth and cellular proliferation as well as modulating host response to infection, injury, and inflammation. Cytokines also influence reproduction and bone remodeling. A large number of cytokines are pleiotropic and share similar functions. In addition, many cytokines influence the production of other cytokines. Analysis and quantification of cytokines in biological fluids has thus become increasingly important. Methods such as bioassay, enzyme-linked immunosorbent assay (ELISA), intracellular staining, ribonuclease protection assay (RPA) and polymerase chain reaction (PCR) have all been used for quantifying cytokines, however, each of these techniques has limitations associated with it. These techniques are not capable of measuring multiple cytokines simultaneously in a limited sample volume.

Assessing the levels of multiple cytokines may be more revealing than analyzing a single protein. Quantifying multiple cytokines on an individual level can be time consuming and expensive. When combined with separately available analyte-specific microparticle sets, this kit is an excellent tool for simultaneously assessing the levels of multiple human cytokines in a single sample. For ease of use, the microparticles are premixed.

Analyte	Microparticle Region
CD40 Ligand	25
EGF	33
Eotaxin	13
FGF basic	77
FIt-3 Ligand	35
Fractalkine	26
G-CSF	36
GM-CSF	37
Granzyme B	38
GRO $\alpha$	27
GRO $\beta$	29
IFN- $\alpha$	39
IFN- $\beta$	42
IFN- $\gamma$	43
IL-1 $\alpha$	78
IL-1 $\beta$	53
IL-1ra	54
IL-2	55
IL-3	56
IL-4	61
IL-5	62
IL-6	63
IL-7	64

Analyte	Microparticle Region
IL-8	30
IL-10	44
IL-12 p70	45
IL-13	46
IL-15	47
IL-17A	48
IL-17E	51
IL-33	57
IP-10	28
MCP-1	15
MIP-1 $\alpha$	20
MIP-1 $\beta$	21
MIP-3 $\alpha$	18
MIP-3 $\beta$	14
PDGF-AA	66
PDGF-AB/BB	67
PD-L1/B7-H1	12
RANTES	72
TGF- $\alpha$	73
TNF- $\alpha$	74
TRAIL	75
VEGF	76

## PRINCIPLE OF THE ASSAY

Luminex® Performance Assay multiplex kits are designed for use with the Luminex® MAGPIX® CCD Imager. Alternatively, kits can be used with the Luminex® 100/200™, Luminex® FLEXMAP 3D®, or Bio-Rad® Bio-Plex®, dual laser, flow-based sorting and detection platforms.

Analyte-specific antibodies are pre-coated onto magnetic microparticles embedded with fluorophores at set ratios for each unique microparticle region. Microparticles, standards and samples are pipetted into wells and the immobilized antibodies bind the analytes of interest. After washing away any unbound substances, a biotinylated antibody cocktail specific to the analytes of interest is added to each well. Following a wash to remove any unbound biotinylated antibody, streptavidin-phycoerythrin conjugate (Streptavidin-PE), which binds to the biotinylated antibody, is added to each well. Final washes remove unbound Streptavidin-PE, the microparticles are resuspended in buffer and read using the Luminex® MAGPIX® Analyzer. A magnet in the analyzer captures and holds the superparamagnetic microparticles in a monolayer. Two spectrally distinct Light Emitting Diodes (LEDs) illuminate the microparticles. One LED excites the dyes inside each microparticle to identify the region and the second LED excites the PE to measure the amount of analyte bound to the microparticle. A sample from each well is imaged with a CCD camera with a set of filters to differentiate excitation levels.

Analysis with the Luminex® 100/200™, Luminex® FLEXMAP 3D®, or Bio-Rad Bio-Plex uses one laser to excite the dyes inside each bead to identify the bead region and the second laser to excite the PE to measure the amount of analyte bound to the bead. All fluorescence emissions from each bead as it passes through the flow cell is then analyzed to differentiate emission levels using a Photomultiplier Tube (PMT) and an Avalanche Photodiode.

## LIMITATIONS OF THE PROCEDURE

- FOR RESEARCH USE ONLY. NOT FOR USE IN DIAGNOSTIC PROCEDURES.
- The kit should not be used beyond the expiration date on the kit label.
- Do not mix or substitute reagents with those from other lots or sources.
- If samples fall outside the dynamic range of the assay, further dilute the samples with the appropriate calibrator diluent and repeat the assay.
- Any variation in diluent, operator, pipetting technique, washing technique, incubation time or temperature, and kit age can cause variation in binding.
- Variations in sample collection, processing, and storage may cause sample value differences.
- This assay is designed to eliminate interference by other factors present in biological samples. Until these factors have been tested in the Luminex® Performance Assay, the possibility of interference cannot be excluded.
- Luminex® Performance Assays afford the user the benefit of multi-analyte analysis of biomarkers in a single complex sample. For each sample type, a single multipurpose diluent is used to optimize recovery, linearity, and reproducibility. Such a multipurpose diluent may not optimize any single analyte to the same degree that a unique diluent selected for analysis of that analyte can optimize conditions, therefore, some performance characteristics may be more variable than those for assays designed specifically for single analyte analysis.
- **Only the analytes listed on the Standard Value Card can be measured with this kit.**

## MATERIALS PROVIDED & STORAGE CONDITIONS

Store the unopened kit at 2-8 °C. Do not use past the kit expiration date.

PART	PART #	DESCRIPTION	STORAGE OF OPENED, DILUTED, OR RECONSTITUTED MATERIAL
Human Performance Panel Standard Cocktail	898855	2 vials of recombinant human cytokines in a buffered protein base with preservatives; 2 lyospheres.	Use a fresh standard and control for each assay. Discard after use.
Human Performance Panel Low Control	898889	2 vials of recombinant human cytokines in a buffered protein base with preservatives; lyophilized.	
Human Performance Panel High Control	898888	2 vials of recombinant human cytokines in a buffered protein base with preservatives; lyophilized.	
Human XL Fixed Panel Microparticle Cocktail	898613	0.600 mL of a concentrated microparticle cocktail with preservative.	May be stored for up to 1 month at 2-8 °C.* <i>Prepare fresh 1X solutions at the time of assay. Discard after use.</i>
Human IL-17E/IL-25 Magnetic Microparticles	898810	0.075 mL a concentrated microparticle stock with preservatives.	
Human CCL5/RANTES Magnetic Microparticles	898840	0.075 mL a concentrated microparticle stock with preservatives.	
Human Performance Panel Biotin-Antibody Cocktail	898854	1 vial of a concentrated biotinylated antibody cocktail; lyophilized.	
Streptavidin-PE	893535	0.250 mL of a concentrated streptavidin-phycoerythrin conjugate with preservatives.	
Microparticle Diluent	895529	6 mL of a buffered protein base with preservative.	May be stored for up to 1 month at 2-8 °C.*
Biotin Antibody Diluent 2	895832	2 vials (5.5 mL/vial) of a buffered protein base with preservative.	
Calibrator Diluent RD6-65	895098	21 mL of a buffered protein base with preservatives. <i>Use diluted 1:2 for cell culture supernate samples. Use undiluted for serum/plasma/platelet-poor plasma samples.</i>	
Wash Buffer Concentrate	895003	21 mL of a 25-fold concentrated solution of buffered surfactant with preservative. <i>May turn yellow over time.</i>	
Microplate	641385	1 flat-bottomed 96-well microplate used as a vessel for the assay.	
Mixing Bottles	895505	2 empty 8 mL bottles used for mixing microparticles with Microparticle Diluent.	
Plate Sealers	640445	4 adhesive foil strips.	
Standard Value Card	750031	1 card listing the standard reconstitution volume and working standard concentrations for this lot of base kit.	
Control Mean Value Card 1	750047	1 card listing the low and high mean control values. <i>For cell culture supernate assay.</i>	
Control Mean Value Card 2	750048	1 card listing the low and high mean control values. <i>For serum/plasma assay.</i>	

\*Provided this is within the expiration date of the kit.

## OTHER SUPPLIES REQUIRED

- **Luminex® Performance Assay analyte-specific kit(s) (see page 1)**
- Luminex® MAGPIX®, Luminex® 100/200™, Luminex® FLEXMAP 3D®, or Bio-Rad Bio-Plex analyzer with X-Y platform
- Hand-held microplate magnet or platewasher with a magnetic platform
- Pipettes and pipette tips
- Deionized or distilled water
- Multi-channel pipette, manifold dispenser, or automated dispensing unit
- 50 mL and 500 mL graduated cylinders
- Horizontal orbital microplate shaker (0.12" orbit) capable of maintaining a speed of  $800 \pm 50$  rpm
- Microcentrifuge
- **Polypropylene** test tubes for dilution of standards and samples

## PRECAUTIONS

Some components in this kit contain a preservative which may cause an allergic skin reaction. Avoid breathing mist.

Wear protective gloves, clothing, eye, and face protection. Wash hands thoroughly after handling. Refer to the SDS on our website prior to use.

## TECHNICAL HINTS

- When mixing or reconstituting protein solutions, always avoid foaming.
- To avoid cross-contamination, change pipette tips between additions of each standard level, between sample additions, and between reagent additions. Also, use separate reservoirs for each reagent.
- To ensure accurate results, proper adhesion of plate sealers during incubation steps is necessary.
- Protect microparticles and Streptavidin-PE from light at all times to prevent photobleaching.

## SAMPLE COLLECTION AND STORAGE

**The sample collection and storage conditions listed below are intended as general guidelines. Sample stability has not been evaluated.**

**Cell Culture Supernates** - Remove particulates by centrifugation and assay immediately or aliquot and store samples at  $\leq -20$  °C. Avoid repeated freeze-thaw cycles.

**Serum** - Use a serum separator tube (SST) and allow samples to clot for 30 minutes at room temperature before centrifuging for 15 minutes at 1000 x g. Remove serum and assay immediately or aliquot and store samples at  $\leq -20$  °C. Avoid repeated freeze-thaw cycles.

**Plasma** - Collect plasma using EDTA or heparin as an anticoagulant. Centrifuge for 15 minutes at 1000 x g within 30 minutes of collection. Assay immediately or aliquot and store samples at  $\leq -20$  °C. Avoid repeated freeze-thaw cycles.

**Platelet-poor Plasma** - Collect plasma on ice using EDTA or heparin as an anticoagulant. Centrifuge at 2-8 °C for 15 minutes at 1000 x g within 30 minutes of collection. For more complete platelet removal, an additional centrifugation step of the separated plasma at 1500 x g for 10 minutes at 2-8 °C is recommended. Assay immediately or aliquot and store samples at  $\leq -20$  °C. Avoid repeated freeze-thaw cycles.

**Note:** Citrate plasma is not validated for use in this assay.

Hemolyzed, icteric, and lipemic samples are not suitable for use in this assay.

Heparin plasma and platelet-poor plasma samples are not suitable for use in the Eotaxin assay.

**Some cytokines may be released upon platelet activation. For example, to measure circulating levels of CD40 Ligand, EGF, GRO $\alpha$ , GRO $\beta$ , PDGF-AA, PDGF-AB/BB, and RANTES, platelet-poor plasma should be used. It should be noted that many protocols for plasma preparation, including procedures recommended by the Clinical and Laboratory Standards Institute (CLSI), result in incomplete removal of platelets or platelet activation. This may cause variable and irreproducible results for assays of factors contained in platelets and released by platelet activation.**

## SAMPLE PREPARATION

**Use polypropylene tubes.**

**Note:** On the day of the assay, ALL fresh and previously frozen serum and plasma samples require centrifugation at 16,000 x g for 4 minutes immediately prior to use or dilution.

Cell culture supernate samples require a 2-fold dilution. A suggested 2-fold dilution is 75  $\mu$ L of sample + 75  $\mu$ L of Calibrator Diluent RD6-65 (diluted 1:2)\*. Mix thoroughly.

Serum, plasma, and platelet-poor plasma samples require a 2-fold dilution. A suggested 2-fold dilution is 75  $\mu$ L of sample + 75  $\mu$ L of Calibrator Diluent RD6-65. Mix thoroughly.

When assaying GRO $\beta$ , PDGF-AA, and PDGF-AB/BB, serum samples must be further diluted 5-fold to a final 10-fold dilution. A suggested 10-fold dilution is 30  $\mu$ L of the 2-fold diluted sample + 120  $\mu$ L of Calibrator Diluent RD6-65. Mix thoroughly.

\*See Reagent Preparation section.

## REAGENT PREPARATION

**Bring all reagents to room temperature before use.**

**Wash Buffer** - If crystals have formed in the concentrate, warm to room temperature and mix gently until the crystals have completely dissolved. Add 20 mL of Wash Buffer Concentrate to 480 mL of deionized or distilled water to prepare 500 mL of Wash Buffer.

**Calibrator Diluent RD6-65 (diluted 1:2) - For cell culture supernate samples only.** Add 10 mL of Calibrator Diluent RD6-65 to 10 mL of deionized or distilled water to prepare 20 mL of Calibrator Diluent RD6-65 (diluted 1:2).


**Low and High Kit Controls - Refer to the vial label for reconstitution volume.** Reconstitute the low and high controls with Calibrator Diluent RD6-65 (diluted 1:2) (*for cell culture supernate samples*) or Calibrator RD6-65 (*for serum/plasma/platelet-poor plasma samples*). Allow the controls to sit for a minimum of 15 minutes with gentle agitation prior to plating.

**Biotin-Antibody Cocktail - Refer to the vial label for reconstitution volume.** Reconstitute the Human Performance Panel Biotin-Antibody Cocktail with Biotin Antibody Diluent 2. Allow the antibody cocktail to sit for a minimum of 20 minutes or a minimum of 5 minutes on a nutator.

**Standard - Refer to the Standard Value Card for the reconstitution volume and assigned values.** Reconstitute the Standard lysospheres with Calibrator Diluent RD6-65 (diluted 1:2) (*for cell culture supernate samples*) or Calibrator RD6-65 (*for serum/plasma/platelet-poor plasma samples*). Allow the standard to sit for a minimum of 20 minutes prior to making dilutions.

**Note:** Do NOT vortex standard cocktail. Gentle agitation should be initiated only after the 20-minute reconstitution step is complete. Standard may have a slightly blue appearance once fully reconstituted.

**Use polypropylene tubes.** Pipette 300  $\mu$ L of the reconstituted standard into the Standard 1 tube. Pipette 200  $\mu$ L of Calibrator Diluent RD6-65 (diluted 1:2) (*for cell culture supernate samples*) or Calibrator RD6-65 (*for serum/plasma/platelet-poor plasma samples*) into the remaining tubes. Use Standard 1 to produce a 3-fold dilution series (below). Refer to the analyte specific datasheets for details. Mix each tube thoroughly before the next transfer. Standard 1 serves as the high standard. The appropriate calibrator diluent serves as the blank.

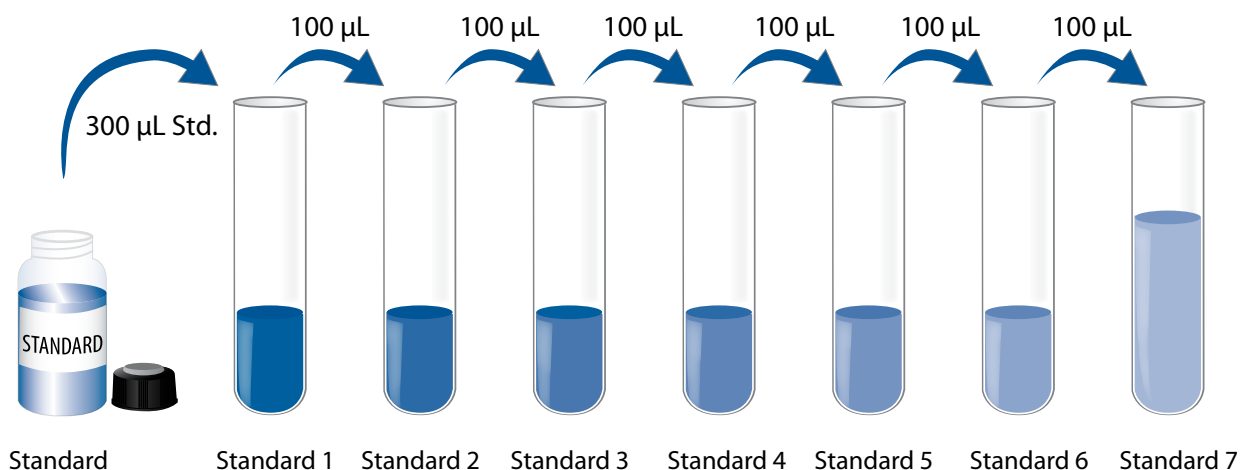



## DILUTED MICROPARTICLE COCKTAIL PREPARATION

1. Dilute the Microparticle Cocktail in the mixing bottle provided.

Number of Wells Used	Microparticle Cocktail	+	Microparticle Diluent
96	0.500 mL	+	5.00 mL
72	0.375 mL	+	3.75 mL
48	0.250 mL	+	2.50 mL
24	0.125 mL	+	1.25 mL

2. To prepare the RANTES and IL-17E microparticles, add the RANTES and IL-17E microparticle stocks into previously diluted microparticles from Step 1 following the table below.

Number of Wells Used	Microparticle Diluent	+	RANTES Microparticles	+	IL-17E Microparticles
96	5.5 mL	+	55 µL	+	55 µL
72	4.125 mL	+	41.25 µL	+	41.25 µL
48	2.75 mL	+	27.5 µL	+	27.5 µL
24	1.375 mL	+	13.75 µL	+	13.75 µL

**Note:** Protect microparticles from light during handling. Diluted microparticles cannot be stored. Prepare microparticles within 30 minutes of use.

## DILUTED BIOTIN-ANTIBODY COCKTAIL PREPARATION

1. Dilute the reconstituted Performance Panel Biotin-Antibody Cocktail in Biotin Antibody Diluent 2. Mix gently.

Number of Wells Used	Biotin-Antibody Cocktail	+	Biotin Antibody Diluent 2
96	0.500 mL	+	5.00 mL
72	0.375 mL	+	3.75 mL
48	0.250 mL	+	2.50 mL
24	0.125 mL	+	1.25 mL

## STREPTAVIDIN-PE PREPARATION

**Use a polypropylene amber bottle or a polypropylene tube wrapped with aluminum foil. Protect Streptavidin-PE from light during handling and storage.**

1. Centrifuge the Streptavidin-PE vial for 30 seconds at 1000 x g prior to removing the cap.
2. Gently vortex the vial, taking precautions not to invert the vial.
3. Dilute the Streptavidin-PE concentrate in Wash Buffer.

Number of Wells Used	Streptavidin-PE Concentrate	+	Wash Buffer
96	220 µL	+	5.35 mL
72	165 µL	+	4.00 mL
48	110 µL	+	2.65 mL
24	55 µL	+	1.35 mL

## INSTRUMENT SETTINGS

**Note:** *Adjust the probe height setting on the analyzer to avoid puncturing the plate. Calibrate the analyzer using the proper reagents for superparamagnetic microparticles (refer to instrument manual).*

### **Luminex® MAGPIX® analyzer:**

- a) Sample volume: 35 µL
- b) Assign the microparticle region for each analyte being measured (see page 1)
- c) 50 count/region
- d) Collect Median Fluorescence Intensity (MFI)

### **Luminex® 100/200™, Luminex® FLEXMAP 3D® and Bio-Rad Bio-Plex analyzers:**

**Note:** *Ensure that the instrument flow rate is set to the default of 60 µL/minute (fast) for all flow based analyzers.*

- a) Sample volume: 50 µL
- b) Bead Type:
  - i. Luminex® 100/200™ and FLEXMAP 3D® select MagPlex
  - ii. Bio-Rad Bio-Plex Manager use Bio-Plex MagPlex Beads (Magnetic)
- c) Doublet Discriminator gates:
  - i. Luminex® 100/200™ and FLEXMAP 3D® set at 8000 and 16,500
  - ii. Bio-Rad Bio-Plex Manager set at 8000 and 23,000
- d) Reporter Gain Setting:
  - i. Luminex® 100/200™ use Default setting
  - ii. Luminex® FLEXMAP 3D® use Standard PMT setting
  - iii. Bio-Rad Bio-Plex Manager use the low RP1 target value for the CAL2 setting
- e) Assign the microparticle region for each analyte being measured (see page 1)
- f) 50 count/region
- g) Collect MFI

## ASSAY PROCEDURE

**Bring all reagents and samples to room temperature before use. It is recommended that all standards, controls, and samples be assayed in duplicate.**

**Note:** *Protect microparticles and Streptavidin-PE from light at all times.*

1. Prepare all reagents, working standards, and samples as directed in the previous sections.
2. Add 50  $\mu$ L of standard, control, or sample\* per well. A plate layout is provided to record standards and samples assayed.
3. Resuspend the diluted Microparticle Cocktail by inversion or vortexing. Add 50  $\mu$ L of the microparticle cocktail to each well of the microplate. Securely cover with a foil plate sealer. Incubate for 2 hours at room temperature on a horizontal orbital microplate shaker (0.12" orbit) set at  $800 \pm 50$  rpm.
4. Using a magnetic device designed to accommodate a microplate, wash by applying the magnet to the bottom of the microplate, allow 1 minute before removing the liquid, filling each well with Wash Buffer (100  $\mu$ L) and allow 1 minute before removing the liquid again. Complete removal of liquid is essential for good performance.  
**Note:** *Do NOT blot; this may cause a loss of microparticles.* Perform the wash procedure three times.  
  
**Note:** *Refer to the magnetic device user manual for proper wash technique using a round bottom microplate.*
5. Add 50  $\mu$ L of diluted Biotin-Antibody Cocktail to all wells. Securely cover with a foil plate sealer and incubate for 1 hour at room temperature on the shaker set at  $800 \pm 50$  rpm.
6. Repeat the wash as in step 4.
7. Add 50  $\mu$ L of diluted Streptavidin-PE to all wells. Securely cover with a foil plate sealer and incubate for 30 minutes at room temperature on the shaker set at  $800 \pm 50$  rpm.
8. Repeat the wash as in step 4.
9. Resuspend the microparticles by adding 100  $\mu$ L of Wash Buffer to each well. Incubate for 2 minutes on the shaker set at  $800 \pm 50$  rpm.
10. Read within 90 minutes using the Luminex® or Bio-Rad analyzer.  
**Note:** *Resuspend microparticles immediately prior to reading by shaking the plate for 2 minutes on the plate shaker set at  $800 \pm 50$  rpm.*

\*Samples require dilution. See Sample Preparation section.

## ASSAY PROCEDURE SUMMARY

**Note:** Protect microparticles and Streptavidin-PE from light at all times.

- ① Prepare all reagents as instructed.  
↓
- ② Add 50  $\mu$ L of standard, control, or sample\* to each well.  
↓
- ③ Add 50  $\mu$ L of diluted Microparticle Cocktail to each well.  
Incubate for 2 hours at RT on a shaker at 800 rpm.  
↓
- ④ Wash by removing the liquid from each well, filling with 100  $\mu$ L Wash Buffer, and removing the liquid again.  
Perform the wash 3 times.  
↓
- ⑤ Add 50  $\mu$ L of diluted Biotin-Antibody Cocktail to each well.  
Cover and incubate for 1 hour at RT on the shaker at 800 rpm.  
↓
- ⑥ Repeat the wash as in step 4.  
↓
- ⑦ Add 50  $\mu$ L of diluted Streptavidin-PE to each well.  
Incubate for 30 minutes at RT on the shaker at 800 rpm.  
↓
- ⑧ Repeat the wash as in step 4.  
↓
- ⑨ Add 100  $\mu$ L of Wash Buffer to each well.  
Incubate for 2 minutes at RT on the shaker at 800 rpm.  
↓
- ⑩ Read within 90 minutes using a Luminex® or Bio-Rad analyzer  
**Note:** Resuspend microparticles immediately prior to reading.

\*Samples require dilution. See Sample Preparation section.

## **CALCULATION OF RESULTS**

Use the Standard concentrations on the Standard Value Card and calculate 3-fold dilutions for the remaining levels. Average the duplicate readings for each standard and sample and subtract the average blank Median Fluorescence Intensity (MFI).

Create a standard curve for each analyte by reducing the data using computer software capable of generating a five parameter logistic (5-PL) curve-fit.

Since samples have been diluted, the concentration read from the standard curve must be multiplied by the dilution factor.

## **CALIBRATION**

This assay is calibrated against highly purified recombinant human cytokines produced at R&D Systems®.

## PRECISION

**Intra-Assay Precision** - is generated from the mean of the %CV's from 40 reportable results across two different concentrations of analytes in a single serum assay.

**Inter-Assay Precision** - is generated from the mean of the %CV's across two different concentrations of analytes across 25 different serum assays.

Analyte	Intra-Assay (%CV)	Inter-Assay (%CV)
B7-H1/PD-L1	1.8	11.3
CCL11/Eotaxin	3.2	10.0
CCL19/MIP-3 $\beta$	2.2	9.7%
CCL2/MCP-1	1.4	8.9
CCL20/MIP-3 $\alpha$	2.4	9.9
CCL3/MIP-1 $\alpha$	1.5	7.3
CCL4/MIP-1 $\beta$	3.0	9.6
CD40 Ligand	3.3	10.6
CX3CL1/Fractalkine	2.7	11.1
CXCL1/GRO $\alpha$	1.0	8.0
CXCL10/IP-10	2.1	8.8
CXCL2/GRO $\beta$	1.2	8.8
CXCL8/IL-8	1.8	9.0
EGF	3.0	9.6
FGF-basic	4.2	12.6
Flt-3 ligand	2.5	10.1
G-CSF	1.8	10.2
GM-CSF	3.1	10.3
Granzyme B	1.8	11.0
IFN- $\alpha$	2.9	9.9
IFN- $\beta$	3.0	11.0
IFN- $\gamma$	2.8	10.4
IL-10	3.3	9.0

Analyte	Intra-Assay (%CV)	Inter-Assay (%CV)
IL-12 p70	2.3	10.5
IL-13	1.6	9.7
IL-15	2.4	9.9
IL-17A	2.5	11.2
IL-17E/IL-25	1.7	9.2
IL-1 $\alpha$	3.0	10.4
IL-1 $\beta$	2.3	10.0
IL-1ra	2.1	10.3
IL-2	2.2	9.2
IL-3	2.5	10.1
IL-33	2.4	12.3
IL-4	3.2	9.7
IL-5	2.2	12.0
IL-6	2.6	9.8
IL-7	2.5	9.8
PDGF-AA	2.8	15.5
PDGF-AB/BB	2.4	12.4
RANTES	4.7	17.4
TGF- $\alpha$	3.0	9.8
TNF- $\alpha$	2.2	8.7
TRAIL	2.0	10.3
VEGF	2.6	10.3

## ACCURACY

**Linearity** - The data represents mean spiked or natural linearity in serum matrix samples. Samples with natural linearity are marked with an asterisk.

Analyte	n	Linearity Average %
B7-H1/PD-L1	4	106
CCL11/Eotaxin	4	92
CCL19/MIP-3 $\beta$	4	123
CCL2/MCP-1*	4	82
CCL20/MIP-3 $\alpha$	4	107
CCL3/MIP-1 $\alpha$	4	77
CCL4/MIP-1 $\beta$	4	103
CD40 Ligand	4	113
CX3CL1/Fractalkine	4	124
CXCL1/GRO $\alpha$	4	80
CXCL10/IP-10	3	110
CXCL2/GRO $\beta$ *	4	114
CXCL8/IL-8	4	118
EGF*	4	110
FGF-basic	4	161
FIt-3 ligand	4	102
G-CSF	4	209
GM-CSF	4	96
Granzyme B	4	99
IFN- $\alpha$	3	98
IFN- $\beta$	4	148
IFN- $\gamma$	4	97
IL-10	4	111

Analyte	n	Linearity Average %
IL-12 p70	4	106
IL-13	4	112
IL-15	3	88
IL-17A	4	110
IL-17E/IL-25	4	122
IL-1 $\alpha$	4	143
IL-1 $\beta$	4	107
IL-1ra*	4	100
IL-2	4	105
IL-3	4	123
IL-33	4	93
IL-4	3	127
IL-5	3	113
IL-6	3	124
IL-7	4	113
PDGF-AA*	4	90
PDGF-AB/BB*	4	96
RANTES*	4	96
TGF- $\alpha$	4	117
TNF- $\alpha$	4	117
TRAIL	3	123
VEGF	4	103

*\*Natural linearity*

## ACCURACY

**Spiked Recovery** - The data represents mean percent recovery of spiked standards ranging from low, medium, and high concentration in serum matrix samples.

Analyte	n	Recovery Average %
B7-H1/PD-L1	4	109
CCL11/Eotaxin	4	107
CCL19/MIP-3 $\beta$	4	95
CCL2/MCP-1	HE	
CCL20/MIP-3 $\alpha$	4	99
CCL3/MIP-1 $\alpha$	4	139
CCL4/MIP-1 $\beta$	4	111
CD40 Ligand	4	100
CX3CL1/Fractalkine	4	88
CXCL1/GRO $\alpha$	HE	
CXCL10/IP-10	HE	
CXCL2/GRO $\beta$	3	106
CXCL8/IL-8	4	96
EGF	4	106
FGF-basic	3	80
Flt-3 ligand	4	109
G-CSF	4	43
GM-CSF	4	110
Granzyme B	4	118
IFN- $\alpha$	4	101
IFN- $\beta$	4	69
IFN- $\gamma$	4	112
IL-10	4	104

Analyte	n	Recovery Average %
IL-12 p70	4	108
IL-13	4	103
IL-15	3	89
IL-17A	4	106
IL-17E/IL-25	4	87
IL-1 $\alpha$	4	80
IL-1 $\beta$	4	99
IL-1ra	4	145
IL-2	4	102
IL-3	4	88
IL-33	4	110
IL-4	3	77
IL-5	3	81
IL-6	3	90
IL-7	4	96
PDGF-AA	HE	
PDGF-AB/BB	HE	
RANTES	HE	
TGF- $\alpha$	4	96
TNF- $\alpha$	4	69
TRAIL	4	88
VEGF	4	109

HE = Spiked recovery was not observed due to natural linearity in high endogenous samples.



## SPECIFICITY

The assay was tested for cross-reactivity and interference with the following factors. Less than 0.5% cross-reactivity and interference was observed unless otherwise noted on the analyte specific datasheet.

### Recombinant human:

4-1BB Ligand	FGF-7/KGF	IL-17A/F	PIGF
6Ckine	FGF-16	IL-17B	SDF-1 $\alpha$
$\alpha_2$ -Macroglobulin	Fibrinogen	IL-17B R	SDF-1 $\beta$
Amphiregulin	G-CSF R	IL-17C	Serpin B9/PI-9
Angiogenin	GITR Ligand	IL-17D	ST2
APRIL	GM-CSF R $\alpha$	IL-17F	Syndecan-2
B7-1	GM-CSF R $\beta$	IL-18	THBS
B7-2	gp130	IL-18 R $\alpha$	TNF RI
B7-H2	GRO $\gamma$	IL-18 R $\beta$	TNF RII
B7-H3	HB-EGF	IL-23p40-p19 het	TRAIL R3
B7-H4	HCC-4	IL-27	TRAIL R4
B7-H6	HGF	IL-36R $\alpha$ /FIL-1 $\delta$	TRANCE
B7-H7	I-309	IL-36 $\alpha$	TSG-6
BAFF/BLyS	IFN- $\alpha$ / $\beta$ R1	IL-36 $\beta$	TGS-14/Pentraxin-3
BLC/BCA-1	IFN- $\alpha$ / $\beta$ R2	IL-36 $\gamma$	TWEAK
Cathepsin C	IFN- $\gamma$ R1	IL-37/FIL-1 $\zeta$ /IL-1F7	VEGF R1/Flt-1
Cathepsin H	IGF-I	IL-38/IL-1F10	VEGF R2/Flk-1/KDR
CCL7/MCP-3	IGF-II	Integrin $\alpha 5\beta 3$	VEGF/PIGF
CCL22/MDC	IL-1 RAcP/IL-1 R3	I-TAC	
CD27 Ligand/CD70	IL-1 RI	LAP (TGF- $\beta$ 1)	
CD30 Ligand	IL-1 RII	LIF	
CNTF	IL-2 R $\alpha$	LIF R	
Cripto-1	IL-3 R $\alpha$	LIGHT	
CXCL5/ENA-78	IL-4 R	LT- $\alpha$ /TNF- $\beta$	
CXCL6/GCP-2/LIX	IL-5 R $\alpha$	LT $\alpha 1/\beta 2$	
CXCL7/NAP-2	IL-6 R $\alpha$	LT $\alpha 2/\beta 1$	
EGFR	IL-6 R $\alpha$ /gp130	MCP-2	
ErbB2	IL-9	MCP-4	
ErbB3	IL-11	M-CSF	
ErbB4	IL-12 R $\beta$ 1	MFG-E8	
Fas Ligand	IL-12 R $\beta$ 2	MIG	
FGF R1 $\alpha$ (IIIb)	IL-12p40	OPG	
FGF R1 $\alpha$ (IIIc)	IL-13 R $\alpha$ 1	OSM	
FGF R $\beta$ (IIIb)	IL-13 R $\alpha$ 2	OX40 Ligand	
FGF R1 $\beta$ (IIIc)	IL-15 R $\alpha$	PARC/MIP-4	
FGF acidic	IL-15/IL-15 R $\alpha$ Complex	PD-1	
FGF-4	IL-17 R	PD-L2	
FGF-5	IL-17 RC	PF-4	
FGF-6	IL-17 RD	Pleiotrophin/PTN	

## SPECIFICITY *CONTINUED*

### Recombinant human multiplex partners:

CCL2/MCP-1	IL-1ra/IL-1F3
CCL3/MIP-1 $\alpha$	IL-2
CCL4/MIP-1 $\beta$	IL-3
CCL5/RANTES	IL-4
CCL11/Eotaxin	IL-5
CCL19/MIP-3 $\beta$	IL-6
CCL20/MIP-3 $\alpha$	IL-7
CD40 Ligand/TNFSF5	IL-8/CXCL8
CXCL1/GRO $\alpha$	IL-10
CXCL2/GRO $\beta$	IL-12 p70
CX3CL1/Fractalkine	IL-13
CXCL10/IP-10	IL-15
EGF	IL-17A
FGF basic	IL-17E/IL-25
Flt-3 Ligand	IL-33
G-CSF	PDGF-AA
GM-CSF	PDGF-AB/BB
Granzyme B	PD-L1/B7-H1
IFN- $\alpha$	TGF- $\alpha$
IFN- $\beta$	TNF- $\alpha$
IFN- $\gamma$	TRAIL/TNFSF10
IL-1 $\alpha$ /IL-1F1	VEGF
IL-1 $\beta$ /IL-1F2	

# PLATE LAYOUT

Use this plate layout to record standards and samples assayed.

12									
11									
10									
9									
8									
7									
6									
5									
4									
3									
2									
1									
	A	B	C	D	E	F	G	H	

**NOTES**

*All trademarks and registered trademarks are the property of their respective owners.*