

# QuantiGlo<sup>®</sup> ELISA

## Human IL-1 $\beta$ /IL-1F2 Immunoassay

Catalog Number QLB00B

For the quantitative determination of human Interleukin 1 beta (IL-1 $\beta$ ) concentrations in cell culture supernates, serum, and plasma.

**Note: The standard reconstitution method has changed. Read this package insert in its entirety before using this product.**

This package insert must be read in its entirety before using this product.  
For research use only. Not for use in diagnostic procedures.

# TABLE OF CONTENTS

SECTION	PAGE
INTRODUCTION .....	1
PRINCIPLE OF THE ASSAY.....	2
TECHNICAL HINTS.....	2
PRECAUTIONS.....	2
MATERIALS PROVIDED & STORAGE CONDITIONS .....	3
OTHER SUPPLIES REQUIRED .....	3
SAMPLE COLLECTION & STORAGE.....	4
SAMPLE PREPARATION.....	4
REAGENT PREPARATION.....	5
ASSAY PROCEDURE .....	6
CALCULATION OF RESULTS.....	7
TYPICAL DATA.....	7
PRECISION .....	8
RECOVERY.....	8
LINEARITY.....	8
SENSITIVITY .....	9
CALIBRATION .....	9
SAMPLE VALUES.....	9
SPECIFICITY.....	10
REFERENCES.....	11
PLATE LAYOUT .....	12

## MANUFACTURED AND DISTRIBUTED BY:

### USA & Canada | R&D Systems, Inc.

614 McKinley Place NE, Minneapolis, MN 55413, USA  
TEL: (800) 343-7475 (612) 379-2956 FAX: (612) 656-4400  
E-MAIL: info@RnDSystems.com

## DISTRIBUTED BY:

### UK & Europe | R&D Systems Europe, Ltd.

19 Barton Lane, Abingdon Science Park, Abingdon OX14 3NB, UK  
TEL: +44 (0)1235 529449 FAX: +44 (0)1235 533420  
E-MAIL: info@RnDSystems.co.uk

### China | R&D Systems China Co., Ltd.

24A1 Hua Min Empire Plaza, 726 West Yan An Road, Shanghai PRC 200050  
TEL: +86 (21) 52380373 FAX: +86 (21) 52371001  
E-MAIL: info@RnDSystemsChina.com.cn

## INTRODUCTION

The Interleukin 1 (IL-1) family of proteins consists of the classic members IL-1 $\alpha$ , IL-1 $\beta$ , and IL-1ra, plus IL-18, IL-33 and IL-1F5-F10. IL-1 $\alpha$  and IL-1 $\beta$  bind to the same cell surface receptors and share biological functions (1). IL-1 is not produced by unstimulated cells of healthy individuals with the exception of skin keratinocytes, some epithelial cells, and certain cells of the central nervous system. In response to inflammatory agents, infections, or microbial endotoxins, however, a dramatic increase in the production of IL-1 by macrophages and various other cell types is observed. IL-1 $\beta$  plays a central role in immune and inflammatory responses, bone remodeling, fever, carbohydrate metabolism, and GH/IGF-I physiology. Inappropriate or prolonged production of IL-1 has been implicated in a variety of pathological conditions including sepsis, rheumatoid arthritis, inflammatory bowel disease, acute and chronic myelogenous leukemia, insulin-dependent diabetes mellitus, atherosclerosis, neuronal injury, and aging-related diseases (2-5).

IL-1 $\alpha$  and IL-1 $\beta$  are structurally related polypeptides that show approximately 25% homology at the amino acid (aa) level. Both are synthesized as 31 kDa precursors that are subsequently cleaved into mature proteins of approximately 17.5 kDa (6, 7). Cleavage of the IL-1 $\beta$  precursor by Caspase-1/ICE is a key step in the inflammatory response (2, 8). Neither IL-1 $\alpha$  nor IL-1 $\beta$  contains a typical hydrophobic signal peptide (9-11), but evidence suggests that these factors can be secreted by non-classical pathways (12, 13). A portion of unprocessed IL-1 $\alpha$  can be presented on the cell membrane and may retain biological activity (14). The precursor form of IL-1 $\beta$ , unlike the IL-1 $\alpha$  precursor, shows little or no biological activity in comparison to the processed form (13, 15). Both unprocessed and mature forms of IL-1 $\beta$  are exported from the cell.

IL-1 $\alpha$  and IL-1 $\beta$  exert their effects through immunoglobulin superfamily receptors that additionally bind IL-1ra. The 80 kDa transmembrane type I receptor (IL-1 RI) is expressed on T cells, fibroblasts, keratinocytes, endothelial cells, synovial lining cells, chondrocytes, and hepatocytes (16, 17). The 68 kDa transmembrane type II receptor (IL-1 RII) is expressed on B cells, neutrophils, and bone marrow cells (18). The two IL-1 receptor types show approximately 28% homology in their extracellular domains but differ significantly in that the type II receptor has a cytoplasmic domain of only 29 aa, whereas the type I receptor has a 213 aa cytoplasmic domain. IL-1 RII does not appear to signal in response to IL-1 and may function as a decoy receptor that attenuates IL-1 function (19). The IL-1 receptor accessory protein (IL-1 RAcP) associates with IL-1 RI and is required for IL-1 RI signal transduction (20). IL-1ra is a secreted molecule that functions as a competitive inhibitor of IL-1 (21, 22). Soluble forms of both IL-1 RI and IL-1 RII have been detected in human plasma, synovial fluids, and the conditioned media of several human cell lines (23, 24). In addition, IL-1 binding proteins that resemble soluble IL-1 RII are encoded by vaccinia and cowpox viruses (25).

The QuantiGlo<sup>®</sup> Human IL-1 $\beta$ /IL-1F2 Chemiluminescent Immunoassay is a 5.5 hour solid phase ELISA designed to measure human IL-1 $\beta$  in cell culture supernates, serum, and plasma. It contains *E. coli*-expressed recombinant human IL-1 $\beta$  and antibodies raised against the recombinant protein. It has been shown to accurately quantitate the recombinant factor. Results obtained using natural IL-1 $\beta$  showed linear curves that were parallel to the standard curves obtained using the QuantiGlo<sup>®</sup> kit standards. These results indicate that this kit can be used to determine relative mass values for natural IL-1 $\beta$ . Reports indicate that this and other ELISA kits calibrated using mature IL-1 $\beta$  as a standard will detect, but considerably underestimate, the unprocessed IL-1 $\beta$  precursor present in samples (26, 27). In biological samples other than cell lysates, the precursor form of IL-1 $\beta$  is usually not the predominant form of IL-1 $\beta$  present and, additionally, is not biologically active. Therefore, results obtained using this kit should provide a useful measure of the levels of active IL-1 $\beta$  present in biological fluids.

## PRINCIPLE OF THE ASSAY

This assay employs the quantitative sandwich enzyme immunoassay technique. A monoclonal antibody specific for human IL-1 $\beta$  has been pre-coated onto a microplate. Standards and samples are pipetted into the wells and any IL-1 $\beta$  present is bound by the immobilized antibody. After washing away any unbound substances, an enzyme-linked polyclonal antibody specific for human IL-1 $\beta$  is added to the wells. Following a wash to remove any unbound antibody-enzyme reagent, an enhanced luminol/peroxide substrate solution is added to the wells and light is produced in proportion to the amount of IL-1 $\beta$  bound in the initial step. A microplate luminometer is used to measure the intensity of the light emitted.

## TECHNICAL HINTS

- FOR RESEARCH USE ONLY. NOT FOR USE IN DIAGNOSTIC PROCEDURES.
- The kit should not be used beyond the expiration date on the kit label.
- Do not mix or substitute reagents with those from other lots or sources.
- If samples generate values higher than the highest standard, dilute the samples with the appropriate calibrator diluent and repeat the assay.
- This assay is designed to eliminate interference by other factors present in biological samples. Until all factors have been tested in the QuantiGlo<sup>®</sup> Immunoassay, the possibility of interference cannot be excluded.
- Variation in pipetting technique, washing technique, luminometers, incubation time or temperature can cause variations in binding.
- Variations in sample collection, processing, and storage may cause sample value differences.
- When mixing or reconstituting protein solutions, always avoid foaming.
- To avoid cross-contamination, change pipette tips between additions of each standard level, between sample additions, and between reagent additions. Also, use separate reservoirs for each reagent.
- To ensure accurate results, proper adhesion of plate sealers during incubation steps is necessary.
- Relative light units (RLUs) may differ among luminometers. Adjust settings as recommended by the instrument manufacturer.
- Relative light units may vary within the 15 minute reading window.
- Due to limitations of many luminometers, it is recommended that the standards be assayed in duplicate from high to low beginning with the high standard in wells A<sub>1</sub> and A<sub>2</sub>.

## PRECAUTIONS

IL-1 $\beta$  is detectable in saliva. Take precautionary measures to prevent contamination of kit reagents while running this assay.

Calibrator Diluent RD6-10 contains sodium azide, which may react with lead and copper plumbing to form explosive metallic azides. Flush with large volumes of water during disposal.

Some components in this kit contain a preservative which may cause an allergic skin reaction. Avoid breathing mist.

Wear protective gloves, clothing, eye, and face protection. Wash hands thoroughly after handling. Refer to the MSDS on our website prior to use.

## MATERIALS PROVIDED & STORAGE CONDITIONS

Store the unopened kit at 2-8 °C. Do not use past kit expiration date.

PART	PART #	DESCRIPTION	STORAGE OF OPENED/ RECONSTITUTED MATERIAL
Human IL-1 $\beta$ Microplate <sup>†</sup>	893171	96 well polystyrene microplate (12 strips of 8 wells) coated with a monoclonal antibody specific for human IL-1 $\beta$ .	Return unused wells to the foil pouch containing the desiccant pack. Reseal along entire edge of the zip-seal. May be stored for up to 1 month at 2-8 °C.*  May be stored for up to 1 month at 2-8 °C.*
Human IL-1 $\beta$ Conjugate <sup>†</sup>	893172	21 mL of a polyclonal antibody specific for human IL-1 $\beta$ conjugated to horseradish peroxidase with preservatives.	
Human IL-1 $\beta$ Standard <sup>†</sup>	890628	Recombinant human IL-1 $\beta$ in a buffered protein base with preservatives; lyophilized. <i>Refer to the vial label for reconstitution volume.</i>	
Assay Diluent RD1W	895117	11 mL of a buffered protein base with preservatives.	
Calibrator Diluent RD6-10	895468	21 mL of buffered protein base with preservatives.	
Wash Buffer Concentrate	895222	100 mL of a 10-fold concentrated solution of buffered surfactant with preservative.	
Glo Reagent A	895868	4 mL of stabilized enhanced luminol.	
Glo Reagent B	895869	8 mL of stabilized hydrogen peroxide.	
Plate Sealers	N/A	4 adhesive strips.	

\* Provided this is within the expiration date of the kit.

<sup>†</sup> This product is covered by the following U.S. patent: 5,681,933.

## OTHER SUPPLIES REQUIRED

- Luminometer set with the following parameters: 1.0 minute lag time; 0.5 sec/well read time; summation mode; auto gain on, or equivalent.
- Pipettes and pipette tips.
- 1000 mL graduated cylinder.
- Deionized or distilled water.
- Squirt bottle, manifold dispenser, or automated microplate washer.
- Horizontal orbital microplate shaker (0.12" orbit) capable of maintaining a speed of 500  $\pm$  50 rpm.
- Test tubes for dilution of standards and samples.
- Human IL-1 $\beta$  Controls (optional; R&D Systems®, QC190).

## SAMPLE COLLECTION & STORAGE

The sample collection and storage conditions listed below are intended as general guidelines. Sample stability has not been evaluated.

**Cell Culture Supernates** - Remove particulates by centrifugation and assay immediately or aliquot and store samples at  $\leq -20$  °C. Avoid repeated freeze-thaw cycles.

**Serum** - Use a serum separator tube (SST) and allow samples to clot for 30 minutes at room temperature before centrifugation for 15 minutes at 1000 x g. Remove serum and assay immediately or aliquot and store samples at  $\leq -20$  °C. Avoid repeated freeze-thaw cycles.

**Plasma** - Collect plasma using EDTA, heparin, or citrate as an anticoagulant. Centrifuge for 15 minutes at 1000 x g within 30 minutes of collection. Assay immediately or aliquot and store samples at  $\leq -20$  °C. Avoid repeated freeze-thaw cycles.

## SAMPLE PREPARATION

Cell culture supernate samples may require dilution to read on the standard curve. If samples are suspected to be  $> 1000$  pg/mL, run both undiluted and diluted in Calibrator Diluent RD6-10. If samples require large dilutions, perform an intermediate dilution with culture media and the final dilution with calibrator diluent.

## REAGENT PREPARATION

**Bring all reagents to room temperature before use.**

**Note:** *IL-1 $\beta$  is detectable in saliva and sweat. Take precautionary measures to prevent contamination of the kit reagents while running the assay.*


**Wash Buffer** - If crystals have formed in the concentrate, warm to room temperature and mix gently until the crystals have completely dissolved. Add 100 mL of Wash Buffer Concentrate to deionized or distilled water to prepare 1000 mL of Wash Buffer.

**Working Glo Reagent** - 1 part Glo Reagent A (4 mL) and 2 parts Glo Reagent B (8 mL) should be mixed together 15 minutes to 4 hours before use in a capped plastic container and protected from light. 100  $\mu$ L of the resultant mixture is required per well.

**Note:** *If running the assay in less than 96 wells, mix appropriate amounts of Glo Reagent A and Glo Reagent B. To assay half a plate (48 wells), mix 2 mL of Glo Reagent A with 4 mL of Glo Reagent B. Working Glo Reagent should be discarded after use.*

**Human IL-1 $\beta$  Standard - Refer to the vial label for reconstitution volume.** Reconstitute the Human IL-1 $\beta$  Standard with deionized or distilled water. This reconstitution produces a stock solution of 25,000 pg/mL. Allow the standard to sit for a minimum of 15 minutes with gentle agitation prior to making dilutions.

Pipette 960  $\mu$ L of Calibrator Diluent RD6-10 into the 1000 pg/mL tube. Pipette 400  $\mu$ L into the remaining tubes. Use the stock solution to produce a 3-fold dilution series (below). Mix each tube thoroughly and change pipette tips between each transfer. The 1,000 pg/mL standard serves as the high standard. Calibrator Diluent RD6-10 serves as the zero standard (0 pg/mL).

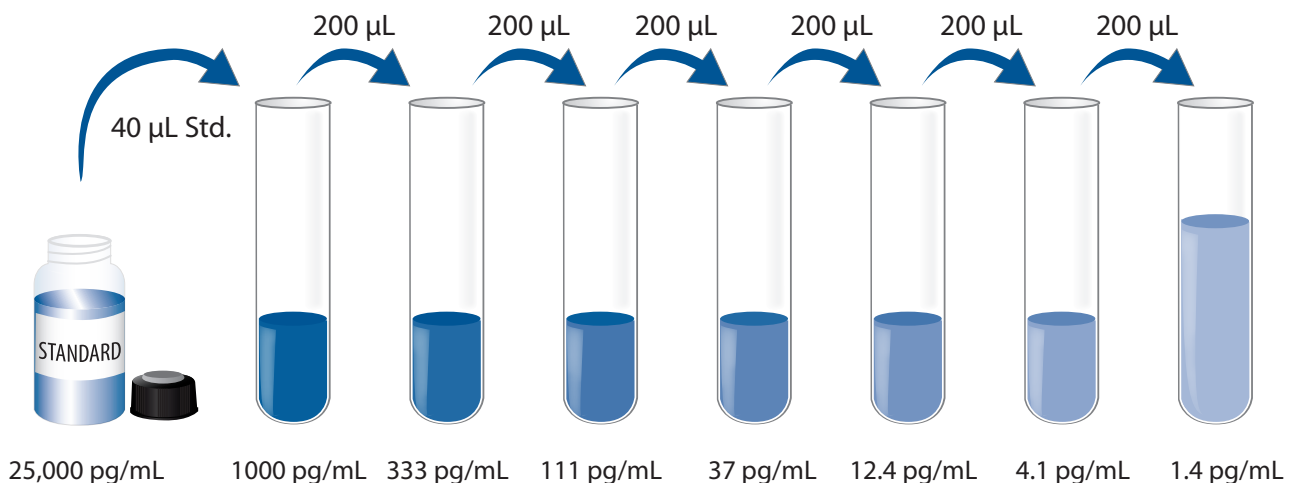

## ASSAY PROCEDURE

**Bring all reagents and samples to room temperature before use. It is recommended that all standards, samples, and controls be assayed in duplicate.**

**Note:** *IL-1 $\beta$  is detectable in saliva and sweat. Take precautionary measures to prevent contamination of the kit reagents while running the assay.*

1. Prepare all reagents, working standards, and samples as directed in the previous sections.
2. Remove excess microplate strips from the plate frame, return them to the foil pouch containing the desiccant pack, and reseal.
3. Add 100  $\mu$ L of Assay Diluent RD1W to each well.
4. Add 100  $\mu$ L of standard, control, or sample per well. Cover with the adhesive strip provided. Incubate for 2 hours at room temperature on a horizontal orbital microplate shaker (0.12" orbit) set at  $500 \pm 50$  rpm. A plate layout is provided for a record of standards and samples assayed.
5. Aspirate each well and wash, repeating the process three times for a total of four washes. Wash by filling each well with Wash Buffer (400  $\mu$ L) using a squirt bottle, manifold dispenser, or autowasher. Complete removal of liquid at each step is essential to good performance. After the last wash, remove any remaining Wash Buffer by aspirating or decanting. Invert the plate and blot it against clean paper towels.
6. Add 200  $\mu$ L of Human IL-1 $\beta$  Conjugate to each well. Cover with a new adhesive strip. Incubate for 3 hours at room temperature on the shaker.  
**Note:** *Prepare Working Glo Reagent at this time.*
7. Repeat the aspiration/wash as in step 5.
8. Add 100  $\mu$ L of Working Glo Reagent to each well. Incubate for 5-20 minutes at room temperature **on the benchtop. Protect from light.**
9. Determine the RLU of each well using a luminometer set with the following parameters: 1.0 min. lag time; 1 sec/well read time; summation mode; auto gain on, or equivalent.

\*Samples may require dilution. See Sample Preparation section.



## CALCULATION OF RESULTS


Average the duplicate readings for each standard, control, and sample and subtract the average zero standard RLU.

Create a standard curve by reducing the data using computer software capable of generating a cubic-spline curve fit.

If samples have been diluted, the concentration read from the standard curve must be multiplied by the dilution factor.

## TYPICAL DATA

This standard curve is provided for demonstration only. A standard curve should be generated for each set of samples assayed.

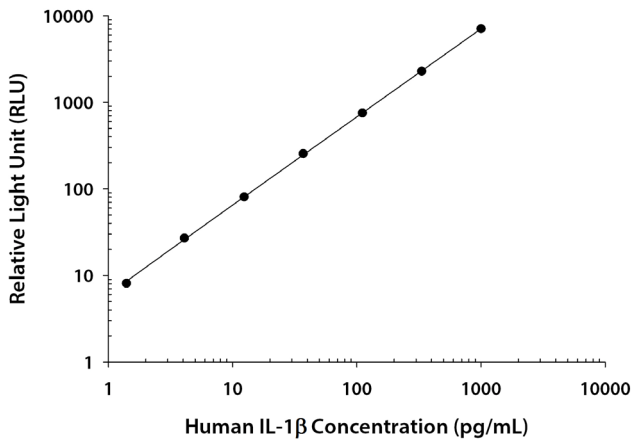

(pg/mL)	RLU	Average	Corrected
0	12.2 12.7	12.4	—
1.4	19.4 21.6	20.5	8.1
4.1	37.0 41.6	39.3	26.9
12.4	91.0 94.0	93.0	80.6
37	261 273	267	255
111	756 770	763	751
333	2292 2298	2295	2283
1000	7085 7085	7085	7073

## PRECISION

### Intra-assay Precision (Precision within an assay)

Four samples of known concentration were tested twenty times on one plate to assess intra-assay precision.

### Inter-assay Precision (Precision between assays)

Four samples of known concentration were tested in twenty separate assays to assess inter-assay precision. Assays were performed by at least three technicians using two lots of components.

Sample	Intra-Assay Precision				Inter-Assay Precision			
	1	2	3	4	1	2	3	4
n	20	20	20	20	20	20	20	20
Mean (pg/mL)	7.96	103	638	4.18	9.86	136	695	5.13
Standard deviation	0.31	5.0	30	0.16	0.71	9.0	38	0.40
CV (%)	3.9	4.9	4.7	3.8	7.2	6.6	5.5	7.8

## RECOVERY

The recovery of recombinant human IL-1 $\beta$  spiked to three different levels in samples throughout the range of the assay in various matrices was evaluated.

Sample Type	Average % Recovery	Range
Cell culture media (n=4)	106	100-115%
Serum (n=4)	96	87-102%
EDTA plasma (n=4)	94	88-107%
Heparin plasma (n=4)	93	87-103%
Citrate plasma (n = 4)	95	90-101%

## LINEARITY

To assess the linearity of the assay, samples containing and/or spiked with high concentrations of human IL-1 $\beta$  in various matrices were diluted with the calibrator diluent to produce samples with values within the dynamic range of the assay.

		Cell culture media (n=4)	Serum (n=4)	EDTA plasma (n=4)	Heparin plasma (n=4)	Citrate plasma (n=4)
1:2	Average % of Expected	98	104	101	101	107
	Range (%)	94-105	102-109	99-104	98-105	103-111
1:4	Average % of Expected	96	105	96	101	111
	Range (%)	92-100	102-108	88-104	95-108	105-116
1:8	Average % of Expected	91	97	95	99	99
	Range (%)	87-96	95-101	87-98	90-106	97-104
1:16	Average % of Expected	97	102	100	103	102
	Range (%)	94-102	97-106	93-104	92-111	100-106
1:32	Average % of Expected	99	102	104	107	104
	Range (%)	92-107	97-107	97-108	95-113	101-104

## SENSITIVITY

Seventy-four assays were evaluated and the minimum detectable dose (MDD) of human IL-1 $\beta$  ranged from 0.05-0.55 pg/mL. The mean MDD was 0.16 pg/mL.

The MDD was determined by adding two standard deviations to the mean RLU of twenty zero standard replicates and calculating the corresponding concentration.

## CALIBRATION

This immunoassay is calibrated against a highly purified *E. coli*-expressed recombinant human IL-1 $\beta$  produced at R&D Systems®.

The NIBSC/WHO IL-1 $\beta$  1st International Standard 86/680, which was intended as a potency standard, was evaluated in this kit. This standard is an *E. coli*-expressed recombinant human IL-1 $\beta$ . The dose response curve of the 1st International Standard parallels the QuantiGlo® standard curve. To convert sample values obtained with the QuantiGlo® Human IL-1 $\beta$  kit to approximate NIBSC 86/680 values, use the equation below.

NIBSC (86/680) approximate value (IU/mL) = 0.0774 x QuantiGlo® Human IL-1 $\beta$  value (pg/mL).

## SAMPLE VALUES

**Serum/Plasma** - Thirty-five serum, EDTA plasma, and heparin plasma sample sets and twenty-one citrate plasma samples drawn from apparently healthy volunteers were evaluated for the presence of human IL-1 $\beta$  in this assay. One serum sample and one EDTA and heparin plasma matched donor sample measured 1.93 pg/mL, 1.99 pg/mL, and 1.99 pg/mL, respectively. All other serum and plasma samples measured below the lowest Human IL-1 $\beta$  Standard. No medical histories were available for the donors used in this study.

**Cell Culture Supernates** - Human peripheral blood mononuclear cells (1 x 10<sup>6</sup> cells/mL) were cultured in RPMI supplemented with 10% fetal bovine serum, 50  $\mu$ M  $\beta$ -mercaptoethanol, and 2 mM L-glutamine. Cells were cultured unstimulated or stimulated with 10  $\mu$ g/mL PHA. Aliquots of the cell culture supernates were removed and assayed for levels of natural human IL-1 $\beta$ .

Condition	Day 1 (pg/mL)	Day 6 (pg/mL)
Stimulated	13.4	3.77
Unstimulated	8120	4940

## SPECIFICITY

This assay recognizes natural and recombinant human IL-1 $\beta$ . This assay also recognizes natural and recombinant rhesus macaque IL-1 $\beta$ .

The factors listed below were prepared at 50 ng/mL in calibrator diluent and assayed for cross-reactivity. Preparations of the following factors at 50 ng/mL in a mid-range recombinant human IL-1 $\beta$  control were assayed for interference. No significant cross-reactivity or interference was observed.

### Recombinant human:

IL-1 $\alpha$   
IL-1F7  
IL-1ra  
IL-1 RAcP  
IL-1 RAPL1  
IL-1 RAPL2  
IL-1 Rrp2  
IL-18  
IL-18 R $\alpha$   
IL-18 R $\beta$   
IL-33  
IL-36 $\alpha$   
IL-36 $\beta$   
IL-36 $\gamma$   
IL-36ra  
IL-38  
SIGIRR  
ST2

### Recombinant mouse:

IL-1 $\alpha$   
IL-1 $\beta$   
IL-33  
SIGIRR  
ST2

### Recombinant rat:

IL-1 $\alpha$   
IL-1ra

### Recombinant porcine:

IL-1 $\alpha$   
IL-1 $\beta$   
IL-1ra

### Recombinant equine:

IL-1 $\beta$   
IL-1ra

### Other recombinants:

cotton rat IL-1 $\alpha$   
canine IL-1 $\beta$   
guinea pig IL-1 $\beta$

Recombinant human Pro-IL-1 $\beta$  cross-reacts approximately 15.4% in this assay.

Recombinant rat IL-1 $\beta$  cross-reacts approximately 1.0% in this assay.

Recombinant cotton rat IL-1 $\beta$  cross-reacts approximately 0.6% in this assay.

## REFERENCES

1. Sims, J.E and D.E. Smith (2010) *Nat. Rev. Immunol.* **10**:89.
2. Martinon, F. and J. Tschopp (2007) *Cell Death Differ.* **14**:10.
3. Isoda, K. and F. Ohsuzu (2006) *J. Atheroscler. Thromb.* **13**:21.
4. Allan, S.M. *et al.* (2005) *Nat. Rev. Immunol.* **5**:629.
5. Kornman, K.S. (2006) *Am. J. Clin. Nutr.* **83**:475S.
6. Giri, J.G. *et al.* (1985) *J. Immunol.* **134**:343.
7. Hazuda, D.J. *et al.* (1990) *J. Biol. Chem.* **265**:6318.
8. Cerretti, D.P. *et al.* (1992) *Science* **256**:97.
9. Lomedico, P.T. *et al.* (1984) *Nature* **312**:458.
10. Auron, P.E. *et al.* (1987) *J. Immunol.* **138**:1447.
11. March, C.J. *et al.* (1985) *Nature* **315**:641.
12. Rubartelli, A. *et al.* (1990) *EMBO J.* **9**:1503.
13. Rubartelli, A. *et al.* (1993) *Cytokine* **5**:117.
14. Kurt-Jones, E.A. *et al.* (1985) *Proc. Natl. Acad. Sci. USA* **82**:1204.
15. Hazuda, D.J. *et al.* (1989) *J. Biol. Chem.* **264**:1689.
16. Urdal, D.L. *et al.* (1988) *J. Biol. Chem.* **263**:2870.
17. Sims, J.E. *et al.* (1988) *Science* **241**:585.
18. McMahan, C.J. *et al.* (1991) *EMBO J.* **10**:2821.
19. Slack, J. *et al.* (1993) *J. Biol. Chem.* **268**:2513.
20. Greenfeder, S.J. *et al.* (1995) *J. Biol. Chem.* **270**:13757.
21. Eisenberg, S.P. *et al.* (1990) *Nature* **343**:341.
22. Carter, D.B. *et al.* (1990) *Nature* **344**:633.
23. Dayer, J-M. and D. Burger (1994) *Eur. Cytokine Netw.* **5**:563.
24. Svenson, M. *et al.* (1993) *Cytokine* **5**:427.
25. Sims, J.E. and S.K. Dower (1994) *Eur. Cytokine Netw.* **5**:539.
26. Herzyk, D.J. *et al.* (1992) *J. Immunol. Methods* **148**:243.
27. Dinarello, C.A. (1992) *J. Immunol. Methods* **148**:255.

## PLATE LAYOUT

Use this plate layout to record standards and samples assayed.

The diagram shows a microplate layout with 12 rows and 8 columns. The rows are numbered 1 through 12 on the left side, and the columns are labeled A through H at the bottom. Each well is represented by a circle. The layout is as follows:

	A	B	C	D	E	F	G	H
1								
2								
3								
4								
5								
6								
7								
8								
9								
10								
11								
12								

**NOTES**

**NOTES**

*All trademarks and registered trademarks are the property of their respective owners.*

©2017 R&D Systems®, Inc.