

## MATERIAL DATA SHEET

### Recombinant Human AQUApure Tetra-Ubiquitin Chains (K63-linked) Cat. # UC-310B

Linkage specific Poly-Ubiquitin chains may be used as a substrate for *in vitro* reactions with deubiquitinating enzymes ("DUB's") that cleave the peptide or isopeptide linkage between adjacent Ubiquitin molecules. Poly-Ubiquitin chains can also be used to investigate mechanisms of binding and recognition between the chains and other proteins that contain Ubiquitin-Associated domains (UBAs), Ubiquitin-interacting motifs (UIMs), ZnF's and/or other Ubiquitin-sensing elements.

K63-linked Tetra-Ubiquitin chains are manufactured using recombinant Ubiquitin and purely enzymatic techniques to avoid the potential for contaminating synthetic intermediates. The correctness of linkage and purity of each production lot is assessed using the Absolute Quantitation of Ubiquitin method (Ub-AQUA), an LCMS-based technique that provides extremely accurate information on the composition of Poly-Ubiquitin samples.

Product Information	
Quantity:	25 μg
MW:	34 kDa
Source:	<i>E. coli</i> -derived human Tetra-Ubiquitin protein Accession # P0CG47
Stock:	1 mg/ml (29 $\mu$ M) in sterile, deionized water
Purity:	>95%, by SDS-PAGE under reducing conditions and visualized by Colloidal Coomassie® Blue stain.
	Ub-AQUA analysis:
	K63: 99.04%
	K11: 0.76%
	K6: 0.13%
	All other linkages $\leq 0.04\%$

Rev. 12/17/2019 Page 1 of 2



# BostonBiochem

An R&D Systems Company

#### **Use & Storage**

Use:	Ubiquitin chains vary in length, linkage, and function. K63-linked Tetra-Ubiquitin Chains (Ub4) are ideal for investigating Ubiquitin-binding proteins and as substrates for Ubiquitin-specific isopeptidases. Reaction conditions will need to be optimized for each specific application. IMPORTANT: Heating this product in SDS-PAGE buffer or terminating reactions containing this product with heated SDS-PAGE buffer could lead to unexpected, high apparent molecular weight banding or smearing on gels that is not representative of product purity. For optimal results, we recommend incubation in SDS-PAGE buffer + DTT at <40 °C for 20 minutes prior to gel electrophoresis.
Storage:	<ul> <li>Use a manual defrost freezer and avoid repeated freeze-thaw cycles.</li> <li>12 months from date of receipt, -20 to -70 °C as supplied.</li> <li>3 months -20 to -70 °C under sterile conditions after opening</li> </ul>

3 months, -20 to -70 °C under sterile conditions after opening.

#### Literature

#### **References:**

- 1. Kirkpatrick D.S. et al. (2006) Nat Cell Biol. 8: 700-10
- 2. Ordureau, A. et al. (2014) Mol. Cell 56: 360-375
- 3. Ordureau, A. et al. (2015) Pro. Nat. Acad. of Sci. USA 112: 6637-6642
- 4. Phu L. et al. (2011) Mol Cell Proteomics 10: M110.003756
- 5. Wall C.E. et al. (2019) Cell Reports 29: 3280

#### For research use only. Not for use in humans.

Rev. 12/17/2019 Page 2 of 2



Global bio-techne.com info@bio-techne.com techsupport@bio-techne.com TEL +1 612 379 2956 USA TEL 800 343 7475 Canada TEL 855 668 8722 China TEL +86 (21) 52380373 Europe | Middle East | Africa TEL +44 (0)1235 529449