

VisUCyte™ HRP Polymer-DAB Cell & Tissue Staining Kit

For the detection of goat, mouse, rabbit, rat, or sheep primary IgG Antibodies with a biotin-free detection system.

Size: 50 Tests

Secondary Antibody-HRP Polymer DAB System

Mouse Kit (Catalog Number VCTS001)

Mouse/Rabbit (Catalog Number VCTS002)

Rabbit Kit (Catalog Number VCTS003)

Goat Kit (Catalog Number VCTS004)

Rat Kit (Catalog Number VCTS005)

Sheep Kit (Catalog Number VCTS006)

**This package insert must be read in its entirety before using this product.
For research use only. Not for use in diagnostic procedures.**

TABLE OF CONTENTS

SECTION	PAGE
INTRODUCTION	1
LIMITATIONS OF THE PROCEDURE	1
TECHNICAL HINTS	2
PRECAUTIONS.....	2
MATERIALS PROVIDED & STORAGE CONDITIONS	3
OTHER SUPPLIES REQUIRED	3
STAINING PROCEDURE.....	4
SUGGESTED READINGS	5

MANUFACTURED AND DISTRIBUTED BY:

USA & Canada | R&D Systems, Inc.

614 McKinley Place NE, Minneapolis, MN 55413, USA
TEL: (800) 343-7475 (612) 379-2956 FAX: (612) 656-4400
E-MAIL: info@RnDSystems.com

DISTRIBUTED BY:

UK & Europe | R&D Systems Europe, Ltd.

19 Barton Lane, Abingdon Science Park, Abingdon OX14 3NB, UK
TEL: +44 (0)1235 529449 FAX: +44 (0)1235 533420
E-MAIL: info@RnDSystems.co.uk

China | R&D Systems China Co., Ltd.

24A1 Hua Min Empire Plaza, 726 West Yan An Road, Shanghai PRC 200050
TEL: +86 (21) 52380373 FAX: +86 (21) 52371001
E-MAIL: info@RnDSystemsChina.com.cn

INTRODUCTION

R&D Systems' VisUCyte™ Cell and Tissue Staining Kits are intended for localization of antigens in a broad range of histological and cytological specimens. These kits may be used with primary IgG antibodies manufactured by R&D Systems or by other vendors. Detection is based on a secondary antibody and HRP enzyme conjugated to a polymer backbone for detection of primary antibodies that react with tissue antigens under study. Visualization is based on enzymatic conversion of a chromogenic substrate 3,3' Diaminobenzidine (DAB) into a colored brown precipitate by horseradish peroxidase (HRP) at the sites of antigen localization, which can then be viewed using bright-field microscopy.

Unlike earlier HRP detection systems, VisUCyte™ kits are biotin-free, eliminating the need to perform biotin and avidin blocking steps in stained samples (eliminates two blocking steps) and using a secondary antibody conjugated to a polymer backbone with a large number of HRP molecules. This one-step detection technique is 3-5 times more sensitive compared to conventional avidin-biotin protocols. VisUCyte™ Polymer detection reagents do not produce non-specific staining which may occur due to binding to phospholipids, nucleic acids and carbohydrate binding proteins.

R&D Systems' VisUCyte™ Cell and Tissue Staining Kits are manufactured to work equally well on cryostat and paraffin-embedded tissue sections, cytopsin preparations (i.e. lymphocytes, monocytes, or transfected cells) and free-floating tissue sections. The kits are made in a ready-to-use format that eliminates the need for extra steps in the procedure, minimizes hands-on time, maximizes convenience and reduces the risk of erroneous calculations.

LIMITATIONS OF THE PROCEDURE

- Do not interchange reagents of this kit with components from other R&D Systems' detection kits and/or similar components manufactured by other vendors.
- Not all primary antibodies are suitable for immunochemistry. Consult your vendor regarding application.
- Any variation in diluents, operator, pipetting technique, washing technique, and incubation time or temperature.
- Avoid dilution of reagents in this kit unless called for in the procedure.
- Reagents should be at room temperature before they are added to samples.

TECHNICAL HINTS

- DAB is not soluble in alcohol or xylene. Xylene-based mounting media may be used to mount tissues stained with DAB.
- Positive Control: To be sure that the staining system is working properly, include a specimen that is known to contain the antigen under study stained by other staining systems (*i.e.*, chromogenic or fluorescent).
- Antibody Positive Control: To be sure that the primary antibodies will detect the antigen under study, include tissues that are known to contain that antigen (*i.e.*, if studying distribution of antigen X in liver that is known to be found in kidney, stain kidney tissue sections along with liver sections).
- Negative Control: This type of control is used to ensure that tissue staining is not caused by non-specific absorption of reagents to the tissue. Three types of negative control may be employed.
 1. **Null Control:** Substitute buffer for primary antibodies. The use of normal serum may not be considered a proper negative control because normal serum may contain unknown antibodies capable of binding to tissue non-specifically that will result in tissue staining.
 2. **Isotype Control:** When using monoclonal primary antibodies incubate control sample with non-immune immunoglobulin of the same isotype.
 3. **Absorption Control:** Incubate with antibodies that were pre-incubated with the corresponding immunogen. Absorption control works satisfactorily if immunogens are peptides. However, if antibodies were raised against proteins, addition of the mixture "antibodies plus proteins" to the tissue may result in higher non-specific staining that is most likely due to the formation of a sandwich complex "tissue-protein-antibody" that is recognized by secondary antibodies. Instead, absorption of antibodies by the immobilized protein may be employed.

PRECAUTIONS

Some components in this kit contain a preservative which may cause an allergic skin reaction. Avoid breathing mist.

3,3' Diaminobenzidine (DAB) is classified as a carcinogen. Wear gloves to avoid DAB contact with skin.

Follow local, state and federal regulations to dispose of used DAB Chromogen.

Wear protective gloves, clothing, eye, and face protection. Wash hands thoroughly after handling. Refer to the MSDS on our website prior to use.

MATERIALS PROVIDED & STORAGE CONDITIONS

Store the unopened kit at 2-8 °C. Do not use past kit expiration date.

PART	PART #	DESCRIPTION	STORAGE OF OPENED/ RECONSTITUTED MATERIAL
Peroxidase Blocking Reagent	865005	6 mL of 3% Hydrogen Peroxide (H ₂ O ₂).	Store at 2-8 °C until expiration date indicated on the kit components.
Mouse IgG Polymer HRP (Vial A)*	898521	6 mL of a anti-mouse secondary antibody in PBS, 1% (w/v) BSA containing a preservative.	
Rabbit IgG Polymer HRP (Vial A)*	898562	6 mL of a anti-rabbit secondary antibody in PBS, 1% (w/v) BSA containing a preservative.	
Mouse/Rabbit IgG Polymer HRP (Vial A)*	898522	6 mL of a anti-mouse/rabbit secondary antibody in PBS, 1% (w/v) BSA containing a preservative.	
Goat IgG Polymer HRP (Vial A)*	898563	6 mL of a anti-goat secondary antibody in PBS, 1% (w/v) BSA containing a preservative.	
Rat IgG Polymer HRP (Vial A)*	898564	6 mL of a anti-rat secondary antibody in PBS, 1% (w/v) BSA containing a preservative.	
Sheep IgG Polymer HRP (Vial A)*	898732	6 mL of a anti-sheep secondary antibody in PBS, 1% (w/v) BSA containing a preservative.	
DAB Chromogen	860001	2 mL of 2.5% 3,3' Diaminobenzidine (DAB) in stabilizing buffer.	
DAB Chromogen Buffer	860005	2 vials (15 mL/vial) of a buffer.	
Serum Blocking Reagent D	865015	6 mL of animal serum in buffer with preservatives.	
Dropper Bottle	720180	15 mL empty dropper bottle.	

*Vial (A) provided is dependent on kit ordered.

OTHER SUPPLIES REQUIRED

- Buffer (PBS, pH 7.4. Tris or Sodium Acetate-based buffers may be substituted).
- Mayer's Hematoxylin Counterstain (Sigma, Catalog # MHS-16) or equivalent.
- Non Aqueous Mounting Medium - Permount™ (Fisher Scientific, Catalog # SP-15-100) or equivalent.

STAINING PROCEDURE

Bring all reagents and samples to room temperature before use. Avoid drying of reagents added to samples by incubating them in a humidity chamber. If drying occurs, discard the specimens and repeat the staining procedure using new samples.

1. Cover sample with 1-3 drops of Peroxidase Blocking Reagent for 5-15 minutes.

Note: *This procedure may affect morphology of unfixed frozen tissue sections. Reduction in incubation time may be required.*

2. Rinse sample with buffer and then wash in buffer for 5 minutes. Wash sample gently since hydrogen peroxide may loosen tissues from the slide.

3. Incubate sample with 1-3 drops of Serum Blocking Reagent D for 15 minutes. Drain slides and carefully wipe off excess Serum Blocking Reagent D before going to the next step.

Do not rinse with buffer.

4. Incubate sample with primary antibody. Follow manufacturer's recommendations regarding working dilution, time and temperature of incubation.

5. Rinse sample with buffer. Wash three times in buffer for 5 minutes/wash.

6. Drain slides and carefully wipe off excess buffer before the next step.

7. Incubate sample with 1-3 drops of Secondary Antibody-polymer HRP (Vial A) for 30-60 minutes. Adjust the incubation time depending on the thickness of the section (the thicker the section, the longer the incubation time) and the affinity of primary antibodies.

8. Repeat step 5.

9. Drain slides and carefully wipe off excess buffer before the next step.

10. Calculate the total volume of DAB Chromogen Solution needed for the entire reaction.

Note: *100-200 μ L of DAB Chromogen Solution is required to cover tissue section on a single slide. Add one drop of DAB Chromogen to 1 mL of DAB Chromogen Buffer.*

11. Mix DAB Chromogen Solution and the DAB Chromogen Buffer in the empty Dropper Bottle provided in the Kit.

12. Add 1-5 drops of freshly prepared DAB Chromogen Solution from the Dropper Bottle to cover the entire sample and incubate for 3-10 minutes. Monitor intensity of staining under a microscope to ensure proper intensity of tissue staining.

Note: *DAB Chromogen Solution is stable for up to six hours.*

13. Rinse with distilled water and then wash in a fresh portion of distilled water for 5 minutes.

Note: *DAB is not soluble in alcohol or xylene. Xylene-based mounting media (i.e. Per-mount) may be used to mount tissues stained with DAB.*

14. Samples stained with DAB may be either mounted without counterstaining or mounted after staining with hematoxylin.

Note: *Excessive counterstaining with hematoxylin may mask areas with weak DAB staining.*

15. Place slides vertically on a filter paper or towel to drain excess mounting medium and let them dry.

16. Slides are ready for observation under the microscope.

SUGGESTED READINGS

1. A.E. Kalyuzhny (2016) *Immunohistochemistry: Essential Elements and Beyond* (Techniques in Life Science and Biomedicine for the Non-Expert); Springer International Publishing AG Part of Springer Nature.
2. Cuello, A.C. ed. (1993) *Immunohistochemistry: Methods in the Neurosciences*, Vol. 14; IBRO Handbook Series, John Wiley & Sons: New York.
3. Bullock, G.R. and P. Petrusz, eds. (1990) *Techniques in Immunocytochemistry*, Vol. 2; Academic Press: New York.
4. Lacey, A.J. ed. (1989) *Light Microscopy in Biology. A Practical Approach*. IRL Press: New York.
5. Bullock, G.R. and P. Petrusz, eds. (1989) *Techniques in Immunocytochemistry*, Vol. 4; Academic Press: New York.
6. Hsu, S.M. *et al.* (1981) *Am. J. Clin. Pathol.* **75**(5):734.
7. Hsu, S.M. *et al.* (1981) *J. Histochem. Cytochem.* **29**(4):577.

All trademarks and registered trademarks are the property of their respective owners.